

electrochemistry handbook



high-quality, user-friendly and cost-effective instruments for every application and budget!



Fisher Scientific puts more into its electrochemical instrument line – so you can get more out of it

As always, Fisher Scientific™ strives to bring you a comprehensive selection of the best laboratory products available. Our updated and expanded selection of electrochemical products features high quality, user-friendly and cost-effective instruments for every application and budget.

The instruments are the state of the art for measuring pH, mV (ORP), ion concentration, conductivity, total dissolved solids (TDS), dissolved oxygen, and temperature. Most of these instruments measure several parameters. The selection guides inside ensure you have the information to make the best choice for your requirements.

Innovative new products, extensive product offering

Fisher Scientific™ accumet™ XL benchtop meters bring advanced features with simplicity to your laboratory. The color touch screen is stunningly vivid and is sure to make a great first impression that doesn't stop. Some newly added features include two stirring probe ports, a 3-position electrode holder, upgradeable software, and USB & RS232 connectivity. We've priced these new meters competitively, so these advanced XL meters have never been more affordable.

Fisher Scientific accumet AB benchtop meters have been the top selling pH meter year after year. That was before we added backlight, multiple views, date/time for GLP requirements, expanded memory, custom buffer calibration options, a calibration alarm, 3-position electrode holder, stirring capability, upgradeable software, and wall-mounting to name a few new features. We have added two new models – the AB250 with ISE capability, and the AB200, our first AB multi-parameter instrument for pH & conductivity. AB basic meters offer great accuracies, simplified performance, and many popular features at an economical price and are poised to continue their success as the top selling pH meter for years to come.

Fisher Scientific accumet waterproof portable meters are the choice of demanding professionals who require rugged, dependable, and feature-rich instruments. These IP67-rated meters deliver exceptional performance and versatility as well as customizable calibration, setup, and measurement capabilities.

Fisher Scientific accumet electrodes meet almost any application. Innovative products like Fisher Scientific™ accuCap™, accu•pHast™ R electrodes, and accuTupH™ XL probes improve performance in tough applications.



AB150
(Page 13)



XL600
(Page 8)



AP125
(Page 18)



13-620-BSP Stirrer
(Page 17)



13-620-632
accuTupH pH/ATC
electrode
(Page 26)



What is Electrochemistry?

The science dealing with the chemical action of electricity and the production of electricity by chemical reactions.

~ wordnetweb.princeton.edu/perl/webwn ~

Benchtop Meters

	Page
Fisher Scientific accumet XL & AB overview and index.....	4
Fisher Scientific accumet XL Series Benchtop Meters.....	6
Fisher Scientific accumet XL600 pH/mV/Temp/ISE/DO/Conductivity Meter.....	8
Fisher Scientific accumet XL500 pH/mV/Temp/ISE/Conductivity Meter.....	10
Fisher Scientific accumet XL200 pH/mV/Temp/Conductivity Meter.....	11
Fisher Scientific accumet XL250 pH/mV/Temp/ISE Meter.....	12
Fisher Scientific accumet XL150 pH/mV/Temp Meter.....	13
Fisher Scientific accumet AB250 pH/mV/ISE/Temp Meter.....	14
Fisher Scientific accumet AB150 pH/mV/Temp Meter.....	15
Fisher Scientific accumet AB200 pH/Conductivity/Temp Meter.....	16
Fisher Scientific accumet Benchtop Stirring Probe.....	17

Portable Meters

Fisher Scientific accumet AP110, AP115 And AP125 Meters.....	18
Fisher Scientific accumet AP71 And AP72 Meters.....	19
Fisher Scientific accumet AP85 And AP75 Meters.....	20
Fisher Scientific accumet AP84 And AP74 Meters.....	21

Electrodes

Fisher Scientific accumet Electrodes.....	22
Fisher Scientific accumet pH Electrodes.....	24
Fisher Scientific accumet pH And Half Cell Electrodes.....	29
Fisher Scientific accumet Electrodes.....	30
Fisher Scientific accumet Half Cell And ORP Electrodes.....	31
Fisher Scientific accumet Ion-Selective Electrodes (ISE).....	32

Probes, And Accessories

Fisher Scientific accumet Conductivity, Dissolved Oxygen And Temperature Probes.....	34
Accessories/Cables/Printers.....	35
Certified Buffer Solutions.....	36

Technical Information

About pH, ORP, And ISE Measurement.....	38
Testing/Care/Troubleshooting.....	40
About Conductivity Measurement.....	42
About Dissolved Oxygen Measurement.....	44

Index.....	46
InnoCal Calibration Services and Certified Fisher Scientific accumet Meters.....	47



The pH Mechanic

– We will keep your lab running smoothly.

Our electrochemistry experts throughout North America will perform a free 18 point inspection of your pH meter and electrodes. Like a trusted mechanic they will diagnose your instruments and recommend parts you need including; calibration settings, electrode check, buffer selection, and more! Subject to availability. To schedule a pH mechanic, contact your Fisher Scientific Sales Representative.

Need more information? Visit www.fishersci.com/accumet for manuals, free communication software, warranty registration, and much more 24/7!



Fisher Scientific accumet High-Quality, Feature-Rich, & User-Friendly Instruments!



Our most popular instruments are now better than ever!

Fisher Scientific strives to bring you a comprehensive selection of the best laboratory products available. Our updated and expanded selection of electrochemistry products offer high quality, user-friendly and cost-effective instruments for every application and budget.

Fisher Scientific accumet was the first to market a touch screen pH meter in 1998 with the AR series, and led the way by introducing the world's first color touch screen pH meter 2003 with the XL series. Our latest Fisher Scientific accumet AB and XL benchtop series once again represent large strides in the world of electrochemistry and stay true to their reputation for having simple-to-understand interfaces.

Whether you are measuring pH, mV (ORP), ion selective concentration, conductivity, total dissolved solids (TDS), dissolved oxygen, temperature, or all of the above, Fisher Scientific accumet has been offering the products you need since 1965. For more information, contact our Fisher Scientific sales representatives around the globe or visit www.fishersci.com.











AB150 Bench Meter
(Page 15)



XL600 Bench Meter
(Pages 8)

Fisher Scientific accumet XL & AB Meter Series

								
Model	XL600	XL500	XL250	XL200	XL150	AB200	AB250	AB150
Page	8	10	12	11	13	16	14	15
Color Touch Screen	✓	✓	✓	✓	✓			
pH-mV (BNC) / Temp	✓✓	✓✓	✓✓	✓	✓	✓	✓	✓
Ion Selective-mV (BNC) / Temp	✓✓	✓✓	✓✓				✓	
Conductivity / TDS / Resistivity / Salinity / Temp	✓	✓		✓		✓		
Dissolved Oxygen / Temp	✓							
Stirring Probe	✓✓	✓✓	✓✓	✓✓	✓	✓	✓	✓
USB Device / RS-232	✓	✓	✓	✓	✓	✓	✓	✓
USB Host / RS-232 / RJ-45 (Ethernet)	✓	✓	✓	✓	✓			

✓✓ = Dual Channel Capacity

Fisher Scientific accumet XL

It's all just a touch away...

Simple Setup Screens

Multi-Channel Measurement

pH Measurement
 Ch 1
7.000
 STABLE
 Last Standardization : 13 Jun 2012 / 2:11:14 AM
 Current Buffer group : USA
 > Touch standardize icon to access standardization mode
 Hide Graph Stop Plotting
 8.00
 7.50
 7.00
 6.50
 0 5 10 15 20 25 30
 Time Elapsed : 0 Hours

Internet Access

Custom Settings

New TFT color touch screen display provides outstanding clarity

Fisher Scientific accumet AB

Better than ever!

New AB backlit graphic display is capable of showing detailed or basic info with a press of a button.

Basic Info

Detailed Info



New Compact Printer
 (Cat. No. 13-637-690) available for direct printing from AB150, AB200, AB250, XL150, XL200, XL250, XL500, or XL600 meters. Includes paper and 100-240 V with 4 plug types for global use.



New 13-637-692 Fisher Scientific Optical USB Mouse for use with any XL series meter. Control your XL meter without touching the screen, or use with your laptop or PC.



New Benchtop Stirring Probe for use with AB or XL series ([Page 35](#))



New 100-240 V Power Supply included with every AB and XL model includes multi-plugs for use around the globe.

Fisher Scientific accumet XL Series Benchtop Meters

Instruments With Features That Will Improve Productivity In Any Laboratory



Removable electrode arm can be mounted on either side of the meter, or add a second arm to facilitate multi-probe operation.



- » TFT color LCD provides crystal clear data from wide viewing angles and at a distance
- » Improved touch screen with glass/glass sensitivity
- » Connect one or two stirring probes; provides mixing without stir plates!
- » Bidirectional communication for LIMS connectivity
- » USB connectivity for use with keyboard, mouse, or other USB device
- » Upgradable application software
- » RJ45 Ethernet/Internet port – connect to Local Area Network (LAN) or Internet, print data to a network printer or surf the web
- » Calibrate with custom buffer values or automatically using one of five preset buffer groups
- » User prompts and context-specific “help” screens guide you through operation
- » Extensive setup screens allow customized settings to meet your specific needs
- » 3-year instrument warranty; register at fishersci.com/accumet
- » Crowded lab bench space? Take advantage of the wall-mounting back panel
- » Do you share your meter with other departments or colleagues? Don't compromise! XL meters allow you to create up to 10 password protected ID's; each includes memory to store 2000 data points - per parameter! Base ID's on individual users, applications, or projects to keep data separate, and retain unique calibration and instrument settings

Choose the minimum or maximum number of channels to display:



Rear View of XL600

STIRRER A:
For use with optional stirring probe (13-620-BSP)

TEMP:
Automatic Temperature Compensation probe (13-620-19) and pH electrodes that also include ATC. *Note: Models XL250, XL500, & XL600 have Dual BNC Channels, each with Temperature*

2/4 CELL:
8-pin DIN connection for 2-cell or 4-cell Cond/TDS/Resistivity/Salinity/Temperature electrodes. *Note: Models XL200, XL500, XL600 only*

POWER:
Power supply, 100/240 VAC, 9 VDC, 18W, center positive. Use only model 13-636-104 (included)


Right



MINI-B USB:
USB communication for field software upgrades (cable included) and data output


RS-232:
RS-232 serial communication output, 2.5 mm jack. For use with RS-232 cable 13-637-681 (included)

Left

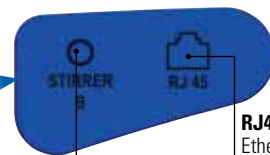

STIRRER B:
For use with optional stirring probe (13-620-BSP)

RJ45:
Ethernet port for network/internet connectivity

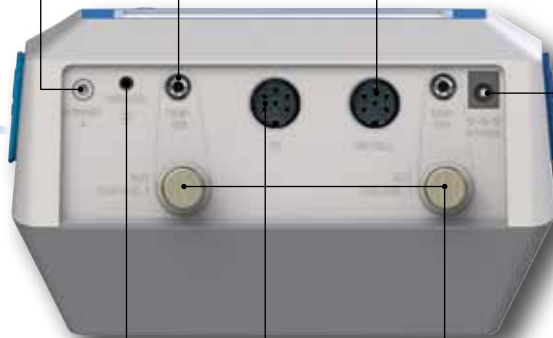

GROUND:
For electrical grounding/electrical noise reduction

DO:
8-pin DIN connection for 13-620-SSP Self-Stirring Dissolved Oxygen/Temperature electrode. *Note: Model XL600 only*

BNC:
For combination pH, mV (ORP, Redox), or Ion Selective Electrodes (ISE). *Note: Models XL250, XL500, & XL600 have Dual BNC Channels, each with Temperature*

Fisher Scientific accumet XL600 pH/mV/Temp/ISE/DO/Conductivity Meter

Quite Possibly The World's Best Meter

- » Customize the XL600 to display 1, 2, 3 or 4 active measurements simultaneously – each with temperature!
Channel 1 : pH, mV, or Ion Selective
Channel 2 : pH, mV, or Ion Selective
Channel 3 : Conductivity, TDS, Resistivity, or Salinity
Channel 4 : Dissolved Oxygen/
BOD/OUR/SOUR
- » Real-time graph provides outstanding detail of any parameter
- » DO mode calculates five-day BOD



*"XL600 Deluxe" kit includes these electrodes.
For ion selective electrodes (ISE), [see page 34](#).*



A complete lab in one instrument. Our top-of-the-line, research-grade touch screen meter measures and displays up to four channels! Unit features two pH/mV/ISE inputs along with one conductivity input (2-cell or 4-cell), as well as one dissolved oxygen input.

Quite simply if you want the best, this is the instrument to put in your lab. Ask your Fisher Scientific representative how you can make your colleagues jealous with accumet, or visit www.fishersci.com/accumet.

For best results, connect optional 13-620-BSP stirrer ([see page 13](#)) for efficient mixing and fast, accurate measurements!

Specifications and Ordering Information

pH	
Range	-2.000 to 20.000
Resolution	0.1/0.01/0.001 selectable
Accuracy	±0.1/0.01/0.002 ±1 LSD
Cal. Points	Up to 6 preset or 5 custom
Buffer Sets	<ul style="list-style-type: none"> • USA: 2.000, 4.010, 6.997, 10.013, 12.000 • NIST: 1.678, 4.010, 6.865, 9.184, 12.460 • DIN: 1.090, 3.060, 4.650, 6.790, 9.230, 12.750 • FSCI: 1.000, 3.000, 6.000, 8.000, 10.000, 13.000 • CUSTOM: Any 2-5 values, ≥1.0 pH unit apart
Slope Display	Yes, with offset
Multiple Slope Display	Yes, up to 5 different slopes
Temp Compensation	Automatic or manual (0 to 100 °C / 32 to 212 °F)
Temp Range (Meter)	0.0 to 100.0 °C / 32.0 to 212.0 °F
Inputs	BNC, ATC
mV	
Range	±2000.0 mV
Rel. mV Range	±2000.0 mV
Resolution	0.1 mV
Accuracy	±0.2 mV or ±0.05 % whichever is greater
Offset Adjustment	Up to ±150 mV
Ion	
Range	0.001 to 19999 (±2000 mV)
Resolution	0.001 / 0.01 / 0.1 / 1 (automatic)
Units	ppm, mg/L, molar
Accuracy	0.5 % full scale (monovalent ion) 1 % full scale (divalent ion)
Cal Points	2 to 6 points from one of following groups; <ul style="list-style-type: none"> • 0.001, 0.01, 0.1, 1, 10, 100 • 0.01, 0.1, 1, 10, 100, 1000 • 0.02, 0.2, 2, 20, 200, 2000 • 0.1, 1, 10, 100, 1000, 10000 • 0.05, 0.5, 5, 50, 500, 5000
Temperature	
Resolution	0.1 °C / 0.1 °F
Accuracy	±0.3 °C / ±0.5 °F
Calibration	Offset in 0.1 ° increments; Offset range: ±5 °C / 9 °F
Conductivity	
Range	0.00 µS to 500.0 mS
Resolution	0.01 / 0.1 µS; 0.001 / 0.01 / 0.1 mS
Accuracy	±1 % full scale
Cal. Points	Automatic (4 points); maximum 1 per range Manual (5 points); maximum 1 per range
Cell Constant	0.010 to 10.000
Cell Types	2 or 4 cell with ATC
Coefficient (Per °C)	Linear & pure; adjustable 0.000 to 10.000 %
Normalization	15.0 to 30.0 °C / 59.0 to 86.0 °F
Compensation	Automatic with supplied cell or manual
Temp Compensation	0.0 to 100 °C / 32.0 to 212.0 °F (0.0 to 80 °C / 32.0 to 176.0 °F with supplied cell)
TDS	
Range	0.00 ppm to 500 ppt (@ TDS factor 1.00)
Resolution	0.01 / 0.1 ppm ; 0.001 / 0.01 / 0.1 ppt
Accuracy	±1 % full scale
Cal. Points	Up to 5
TDS Factor	0.400 to 1.000

Salinity	
Range	0 to 80.0 ppt
Resolution	0.01 / 0.1 ppm; 0.001 / 0.01 / 0.1 ppt
Accuracy	±1 % full scale
Resistivity	
Range	2.000 Ω to 20.0 MΩ
Resolution	0.01 / 0.1 Ω; 0.001 / 0.1 kΩ; 0.01 MΩ
Accuracy	±1 % full scale
Other	
Display	Color TFT Touchscreen, 130 x 100mm (active area)
Output	RJ45, RS232, USB-A, mini-B USB, stirrer port (2)
Adjustable Stirrer Speed	Yes. Each XL can accept (2) stirrers with individual speed control
Language selection	English, French, Spanish, German, Italian, Chinese, Korean, & Portuguese
Memory	2,000 data points per parameter, per user ID, up to 10 user ID's + Default
Datalogging	Manual, timed selectable every 3 to 86400 sec. (24 hours)
Cal Due Alarm	Yes
High / Low Alarms	Yes, user selectable, visual & audible, parameter based
Power	9 V DC adapter, center positive, 1.3 A (100 / 240 VAC, SMPS)

XL600 Deluxe Kit: Includes meter, self-stirring BOD probe (13-620-SSP), TRIS compatible accuTupH pH electrode (13-620-183A), ATC probe (13-620-19), Conductivity/Temp probe (13-620-100), electrode arm, RS-232 & USB cables, 100/240 V power supply, and manual.

XL600 Kit: Includes meter, TRIS compatible accuTupH pH electrode (13-620-183A), ATC probe (13-620-19), Conductivity/Temp probe (13-620-100), electrode arm, RS-232 & USB cables, 100/240 V power supply, and manual.

XL600 Meter Only: Includes meter, electrode arm, RS-232 & USB cables, 100/240 V power supply, and manual.

Description	Catalog No.
XL600 Deluxe Kit	13-636-XL600D
XL600 Kit	13-636-XL600
XL600 Meter Only	13-636-XL600A
Fisher Scientific accumet Benchtop Stirring Probe	13-620-BSP

Fisher Scientific accumet XL500 pH/mV/Temp/ISE/Conductivity Meter

Multi-Channel Operation Simplifies Your Lab Work

- » Up to three channels measured and displayed simultaneously
- » TFT color LCD provides crystal clear data from wide viewing angles and at a distance
- » Improved touch screen with glass/glass sensitivity
- » Connect one or two stirring probes; provides mixing without stir plates!
- » Bidirectional communication for LIMS connectivity
- » USB connectivity for use with keyboard, mouse, or other USB device
- » Upgradable application software
- » RJ45 Ethernet/Internet port – connect to Local Area Network (LAN) or Internet, print data to a network printer or surf the web
- » Calibrate with custom buffer values or automatically using one of five preset buffer groups



Full-function pH and mV capabilities. Extended pH range, with resolution adjustable to 0.001 and calibration from 5 preset buffer groups, custom buffers, and up to six points for exceptional ± 0.002 accuracy.

Ion-selective electrode measurement is easy, with the ion name displayed right on the measurement screen so you'll never have the wrong ISE curve. Automatically run direct or addition/subtraction methods. On-screen prompts allow easy entry of volumes and concentrations,

quickly providing sample concentration without tedious paper calculations.

Measure conductivity, salinity, or resistivity from 0 to 500 mS. Selectable temperature coefficients, cell constants, and display units ensure accurate readings throughout a broad range of samples.

Specifications and Ordering Information

pH Mode	
Range	-2.000 to 20.000
Resolution	0.1 / 0.01 / 0.001
Accuracy	$\pm 0.1 / 0.01 / 0.002 \pm 1$ LSD
Auto Buffer Recognition	5 preset buffer groups + custom
mV Mode	
Range	± 2000 mV / Rel mV
Resolution	0.1
Accuracy	± 0.2
Temperature Mode	
Range	-5.0 to 105.0 °C
Resolution	0.1 °C (0.1 °F)
Accuracy	± 0.2 °C (± 0.3 °F)
Ion Mode	
Range	1×10^{-6} to 9.99×10^{10}
Resolution	0.1 / 0.01 / 0.001
Accuracy	± 0.5 % full scale (monovalent ion) 1 % full scale (divalent ion)
Conductivity Mode	
Range	0 to 500.0 mS
Resolution	0.01 / 0.1 μ S ; 0.001 / 0.01 / 0.1 mS
Accuracy	± 1 % full scale
TDS Mode	
Range	0.050 ppm to 500 ppt (@ TDS factor 1.00)
Salinity Mode	
Range	0 to 80.0 ppt
Resistivity Mode	
Range	2.000 Ω to 20.0 M Ω
Output	RJ45, RS232, USB-A, mini-B USB, stirrer port (2)
Language Selection	English, French, Spanish, German, Italian, Chinese, Korean, & Portuguese
High / Low Alarms	Yes, user selectable, visual & audible, parameter based

XL500 Kit: Includes meter, TRIS compatible accuTupH pH electrode (13-620-183A), ATC probe (13-620-19), Conductivity/Temp probe (13-620-100), electrode arm, RS-232 & USB cables, 100/240 V power supply, and manual.

XL500 Meter Only: Includes meter, electrode arm, RS-232 & USB cables, 100/240 V power supply, and manual.

Description	Catalog No.
XL500 Kit	13-636-XL500
XL500 Meter Only	13-636-XL500A
Fisher Scientific accumet Benchtop Stirring Probe	13-620-BSP



Kit includes 13-620-100 cell.
For other conductivity cells, see page 34.

Fisher Scientific accumet XL200 pH/mV/Temp/Conductivity Meter



No Need To Swap Probes – Measure And Display Two Channels Simultaneously

- » Measure pH/Temp and Conductivity/Temp simultaneously
- » TFT color LCD provides crystal clear data from wide viewing angles and at a distance
- » Improved touch screen with glass/glass sensitivity
- » Connect one or two stirring probes; provides mixing without stir plates!
- » Bidirectional communication for LIMS connectivity
- » USB connectivity for use with keyboard, mouse, or other USB device
- » Upgradable application software
- » RJ45 Ethernet/Internet port – connect to Local Area Network (LAN) or Internet, print data to a network printer or surf the web
- » Calibrate with custom buffer values or automatically using one of five preset buffer groups



Specifications and Ordering Information

pH Mode	
Range	-2.000 to 20.000
Resolution	0.1 / 0.01 / 0.001
Accuracy	±0.1 / 0.01 / 0.002 ±1 LSD
Auto Buffer Recognition	5 preset buffer groups + custom
mV Mode	
Range	±2000 mV / Rel mV
Resolution	0.1
Accuracy	±0.2
Temperature Mode	
Range	-5.0 to 105.0 °C
Resolution	0.1 °C (0.1 °F)
Accuracy	±0.2 °C (±0.3 °F)
Conductivity Mode	
Range	0 to 500.0 mS
Resolution	0.01 / 0.1 µS ; 0.001 / 0.01 / 0.1 mS
Accuracy	±1 % full scale
TDS Mode	
Range	0.050 ppm to 500 ppt (@ TDS factor 1.00)
Salinity Mode	
Range	0 to 80.0 ppt
Resistivity Mode	
Range	2.000 Ω to 20.0 MΩ
Output	RJ45, RS232, USB-A, mini-B USB, stirrer port (2)
Language Selection	English, French, Spanish, German, Italian, Chinese, Korean, & Portuguese
High / Low Alarms	Yes, user selectable, visual & audible, parameter based

XL200 Kit: Includes meter, TRIS compatible accuTupH pH electrode (13-620-183A), ATC probe (13-620-19), Conductivity/Temp probe (13-620-100), electrode arm, RS-232 & USB cables, 100/240 V power supply, and manual.

XL200 Meter Only: Includes meter, electrode arm, RS-232 & USB cables, 100/240 V power supply, and manual.

Description	Catalog No.
XL200 Kit	13-636-XL200
XL200 Meter Only	13-636-XL200A
Fisher Scientific accumet Benchtop Stirring Probe	13-620-BSP



Kit includes 13-620-100 cell.
For other conductivity cells, [see page 34](#).

The XL200 combines two instruments in one; a pH/mV/Temperature meter and Conductivity/TDS/Resistivity/Salinity meter. Display one parameter for better focus and live graphing, or choose dual display with a touch of the screen.

Use one or two optional 13-620-BSP stirring probes for best results. Use optional 13-637-671 to add a second electrode arm and measure on both sides of the XL200.

The XL200 is compatible with a variety of Fisher Scientific accumet conductivity cells to match your requirements. Selectable cell constants let you use K = 1 cells for general purpose, 0.1 cells for high purity water, or 10 cells for brines or other concentrated solutions.



Fisher Scientific accumet XL250 pH/mV/Temp/ISE Meter

Multiple BNC Ports Allow Flexibility

- » Dual-BNC channels; use them for pH or mV or ion selective measurement – you decide
- » Accepts any pH, ORP, or ion selective electrode with BNC connector
- » View independent electrode slope in each channel and each range at a glance
- » TFT color LCD provides crystal clear data from wide viewing angles and at a distance
- » Connect one or two stirring probes; provides mixing without stir plates!
- » Bidirectional communication for LIMS connectivity
- » USB connectivity for use with keyboard, mouse, or other USB device
- » RJ45 Ethernet/Internet port – connect to Local Area Network (LAN) or Internet, print data to a network printer or surf the web
- » Calibrate with custom buffer values or automatically using one of five preset buffer groups

For best results, connect optional 13-620-BSP stirrer (see page 13) for efficient mixing and fast, accurate measurements!

Ion-selective electrode measurement has never been this easy. And with dual-channel capability, there is no need to recalibrate every time you switch from pH to ISE measurement. The XL250 displays the ion name right on the measurement screen so you'll never have the wrong ISE curve.

Addition/subtraction methods are almost foolproof. The XL250 prompts you through the method steps, allowing easy entry of volumes and concentrations, providing you with the sample concentration without tedious paper calculations.

Specifications and Ordering Information

pH Mode	
Range	-2.000 to 20.000
Resolution	0.1 / 0.01 / 0.001
Accuracy	±0.1 / 0.01 / 0.002 ±1 LSD
Auto Buffer Recognition	5 preset buffer groups + custom
mV Mode	
Range	±2000 mV / Rel mV
Resolution	0.1
Accuracy	±0.2
Temperature Mode	
Range	-5.0 to 105.0 °C
Resolution	0.1 °C (0.1 °F)
Accuracy	±0.2 °C (±0.3 °F)
Ion Mode	
Range	1 x 10 ⁻⁶ to 9.99 x 10 ¹⁰
Resolution	0.1 / 0.01 / 0.001
Accuracy	±0.5 % full scale (monovalent ion) 1 % full scale (divalent ion)
Output	RJ45, RS232, USB-A, mini-B USB, stirrer port (2)
Language Selection	English, French, Spanish, German, Italian, Chinese, Korean, & Portuguese
High / Low Alarms	Yes, user selectable, visual & audible, parameter based

Fluoride Kit: Includes meter, 13-620-629 fluoride electrode, TRIS compatible accuTupH pH electrode (13-620-183A), ATC probe (13-620-19), electrode arm, RS-232 & USB cables, 110/220 V power supply, and manual.

Ammonia Kit: Includes meter, 13-620-509 ammonia electrode, TRIS compatible accuTupH pH electrode (13-620-183A), ATC probe (13-620-19), electrode arm, RS-232 & USB cables, 110/220 V power supply, and manual.

Meter Kit: Includes meter, TRIS compatible accuTupH pH electrode (13-620-183A), ATC probe (13-620-19), electrode arm, RS-232 & USB cables, 110/220 V power supply, and manual.

Meter Only: Includes meter, electrode arm, RS-232 & USB cables, 100/240 V power supply, and manual.

Description	Catalog No.
XL250 Fluoride Kit	13-636-XL250F
XL250 Ammonia Kit	13-636-XL250N
XL250 Kit	13-636-XL250
XL250 Meter Only	13-636-XL250A

Ion Selective Electrodes (ISE)

Description	Catalog No.
Fluoride	13-620-629
Ammonia	13-620-509
Nitrate	13-620-534
Chloride	13-620-627
Sodium	13-620-503A



Fisher Scientific accumet XL150 pH/mV/Temp Meter



Simple Yet Powerful, Large Full-Color Touch Screen For Easy Operation

- » Calibrate with custom pH buffer values that you choose or automatically from one of six preset buffer groups including a new pure water buffer group
- » New high-resolution color display provides crystal clear data from wide viewing angles and at a distance
- » Connect one or two stirring probes; provides mixing without stir plates!
- » New bidirectional communication for LIMS connectivity and 3rd party software
- » USB host connectivity for peripherals such as keyboard, mouse, or other USB device
- » RJ45 Ethernet/Internet port – connect to Local Area Network (LAN) or Internet, print data to a network printer or surf the web
- » Optional calibration reminder & high/low alarm
- » Create up to 10 password protected ID's; each includes memory to store 2000 data points. Base ID's on individual users, applications, or projects to keep data separate, and retain unique calibration and instrument settings
- » 3-year instrument warranty; register at fishersci.com/accumet

Specifications and Ordering Information

pH Mode	
Range	-2.000 to 20.000
Resolution	0.1 / 0.01 / 0.001
Accuracy	±0.1 / 0.01 / 0.002 ±1 LSD
Auto Buffer Recognition	USA: 2.00, 4.01, 7.00, 10.01, 12.00 NIST: 1.68, 4.01, 6.86, 9.18, 12.46 DIN 19267: 1.09, 3.06, 4.65, 6.79, 9.23, 12.75 FSCI: 1.00, 3.00, 6.00, 8.00, 10.00, 13.00 Pure Water: 4.10, 6.97, 9.15 CUSTOM: 2-5 points; any values ≥1.0 pH units apart
mV Mode	
Range	±2000 mV / Rel mV
Resolution	0.1
Accuracy	±0.2
Temperature Mode	
Range	-5.0 to 105.0 °C
Resolution	0.1 °C (0.1 °F)
Accuracy	±0.2 °C (±0.3 °F)
Output	RJ45, RS232, USB-A, mini-B USB, stirrer port (2)
Language Selection	English, French, Spanish, German, Italian, Chinese, Korean, & Portuguese
High / Low Alarms	Yes, user selectable, visual & audible, parameter based

XL150 Kit: Includes meter, TRIS compatible accuTupH pH electrode (13-620-183A), ATC probe (13-620-19), electrode arm, RS-232 & USB cables, 110/220 V power supply, and manual.

XL150 Meter Only: Includes meter, electrode arm, RS-232 & USB cables, 100/240 V power supply, and manual.

Description	Catalog No.
XL150 Kit	13-636-XL150
XL150 Meter Only	13-636-XL150A
Fisher Scientific accumet Benchtop Stirring Probe	13-620-BSP

For best results, connect optional 13-620-BSP stirrer (see page 13) for efficient mixing and fast, accurate measurements!



Measure pH and mV with all the power of the top-of-the-line XL600. The XL150 offers simplified, single-channel operation but doesn't skimp on performance or features. Offers the same extended pH range (-2 to 20), with resolution

adjustable to 0.001 and calibration at up to six points for exceptional accuracy (±0.002).

Extensive GLP capabilities with customizable printer or computer output and 2000-reading

datalog capability. Print setup allows you to print exactly what you want, exactly when you want to.



Fisher Scientific accumet AB250 pH/mV/ISE/Temp Meter

Versatile BNC Connection For pH, Redox/ORP, Or Ion Selective Measurements

- » NEW model for pH/mV/Ion Selective measurements
- » Date and Time to meet GLP
- » Password protection
- » Calibration reminder & high/low alarm
- » Toggle views with a touch of a button
- » Graphic LCD with adjustable backlight
- » Calibrate with custom pH buffer values or preset buffer groups
- » 500 point memory – log manually or at intervals
- » Easily view calibration data
- » Upgradable software
- » Selectable 0.1 / 0.01 / 0.001 pH resolution



Affordable Ion measurements have never been easier with the AB250. Connect a pH electrode for use as a pH meter or connect any Ion Selective Electrode when ion measurement is desired.

View detailed calibration data with touch of a button. Collapse the display to show the main reading with temperature, or expand to display additional details.

Specifications and Ordering Information

pH Mode	
Range	-2.000 to 20.000 pH
Resolution	0.1 / 0.01 / 0.001 pH
Accuracy	±0.002 pH + 1 LSD
Buffer Sets	USA: 2.000, 4.010, 6.997, 10.013, 12.000 NIST: 1.678, 4.010, 6.865, 9.184, 12.460 DIN: 1.090, 3.060, 4.650, 6.790, 9.230, 12.750 FSCI: 1.000, 3.000, 6.000, 8.000, 10.000, 13.000 CUSTOM: Any 2-5 values, ≥1.0 pH unit apart
Slope Display	Yes, up to 5 different slopes with offset
Temp Compensation	Automatic or manual (0 to 100 °C / 32 to 212 °F)
mV Mode	
Range	±2000.0 mV / Rel. mV
Resolution	0.1 mV
Accuracy	±0.2 mV or ±0.05 % whichever is greater
Offset Adjustment	Up to ±150 mV
Ion Mode	
Range	0.001 to 19999 (±2000 mV)
Resolution	0.001 / 0.01 / 0.1 / 1 (automatic)
Units	ppm, mg/L, molar
Accuracy	0.5 % full scale (monovalent ion) 1 % full scale (divalent ion)
Cal Points	2 to 6 points from one of following groups; 0.001, 0.01, 0.1, 1, 10, 100 0.01, 0.1, 1, 10, 100, 1000 0.02, 0.2, 2, 20, 200, 2000 0.1, 1, 10, 100, 1000, 10000 0.05, 0.5, 5, 50, 500, 5000
Temperature Mode	
Resolution	0.1 °C / 0.1 °F
Accuracy	±0.3 °C / ±0.5 °F
Calibration	Offset in 0.1 ° increments; offset range: ±5 °C / 9 °F
Output	RS-232 (phono plug), mini-B USB, stirrer
Language Selection	English, French, Spanish, German, Italian, Chinese, Korean, & Portuguese
Memory	500 data sets, viewable
Datalogging	Manual, timed (selectable every 3 to 3600 seconds) Printer or CSV format
Cal Due Alarm	Yes, user selectable from 8 hrs, 16 hrs, & 1-31 days
High / Low Alarms	Yes, user selectable, visual & audible, parameter based

AB250 Kit: Includes meter, TRIS compatible pH/ATC electrode (13-620-631), electrode arm, RS-232 & USB cables, 100/240 V power supply, and manual.

AB250 Meter Only: Includes meter, electrode arm, RS-232 & USB cables, 100/240 V power supply, and manual.

Description	Catalog No.
AB250 Kit	13-636-AB250
AB250 Meter Only	13-636-AB250A

Ion Selective Electrodes (ISE)

Description	Catalog No.
Fluoride	13-620-629
Ammonia	13-620-509
Nitrate	13-620-534
Chloride	13-620-627
Sodium	13-620-503A

For best results, connect optional 13-620-BSP stirrer (see page 13) for efficient mixing and fast, accurate measurements!

Fisher Scientific accumet AB150 pH/mV/Temp Meter

Intuitive, Simple Operation And High Accuracy In A Compact, Affordable Meter



- » Easy-to-understand prompts
- » New graphic display with adjustable backlight
- » Calibrate with custom pH buffer values
- » Date and Time to meet GLP
- » Password protection
- » Optional calibration reminder & high/low alarm
- » 500 point memory – log manually or at intervals
- » Easily view calibration data
- » 0.001 pH resolution and ± 0.002 pH accuracy!



Specifications and Ordering Information

pH Mode	
Range	-2.000 to 20.000 pH
Resolution	0.1 / 0.01 / 0.001 pH
Accuracy	± 0.002 pH + 1 LSD
Buffer Sets	USA: 2.000, 4.010, 6.997, 10.013, 12.000 NIST: 1.678, 4.010, 6.865, 9.184, 12.460 DIN: 1.090, 3.060, 4.650, 6.790, 9.230, 12.750 FSCI: 1.000, 3.000, 6.000, 8.000, 10.000, 13.000 CUSTOM: Any 2-5 values, ≥ 1.0 pH unit apart
Slope Display	Yes, up to 5 different slopes with offset
Temp Compensation	Automatic or manual (0 to 100 °C / 32 to 212 °F)
mV Mode	
Range	± 2000.0 mV / Rel. mV
Resolution	0.1 mV
Accuracy	± 0.2 mV or ± 0.05 % whichever is greater
Offset Adjustment	Up to ± 150 mV
Temperature Mode	
Resolution	0.1 °C / 0.1 °F
Accuracy	± 0.3 °C / ± 0.5 °F
Calibration	Offset in 0.1 ° increments; offset range: ± 5 °C / 9 °F
Output	RS-232 (phono plug), mini-B USB, stirrer
Language Selection	English, French, Spanish, German, Italian, Chinese, Korean, & Portuguese
Memory	500 data sets, viewable
Datalogging	Manual, timed (selectable every 3 to 3600 seconds) Printer or CSV format
Cal Due Alarm	Yes, user selectable from 8 hrs, 16 hrs, & 1-31 days
High / Low Alarms	Yes, user selectable, visual & audible, parameter based

AB150 BioBasic Kit: Includes meter, TRIS compatible accuTupH pH electrode (13-620-183A), ATC probe (13-620-19), electrode arm, 110/220 V power supply, and manual.

AB150 Kit: Includes meter, TRIS compatible pH/ATC electrode (13-620-631), electrode arm, RS-232 & USB cables, 100/240 V power supply, and manual.

AB150 Meter Only: Includes meter, electrode arm, RS-232 & USB cables, 100/240 V power supply, and manual.

Description	Catalog No.
AB150 BioBasic Kit	13-636-AB150B
AB150 Kit	13-636-AB150
AB150 Meter Only	13-636-AB150A

Replacement Parts And Accessories For AB Series Benchtop Meters

Description	Catalog No.
Fisher Scientific accumet™ Benchtop Stirring Probe	13-620-BSP
pH Electrode, Double Junction, Glass Body, Refillable	13-620-183A
pH/ATC Electrode, Double Junction, Plastic Body, Refillable	13-620-631
Temperature Probe, Stainless Steel	13-620-19
Compact Printer, 100-240 V. Includes 1 Roll of Paper	13-637-690
Replacement Power Supply, 100/240 V	13-636-104

For laboratories that need a basic, accurate, easy-to-use pH/mV meter with exceptionally high accuracy and resolution, the AB150 meter is the perfect choice.

Easy operation, accuracy, and affordability set the Fisher Scientific accumet AB150 meter apart from the competition.

Kits include electrodes suitable for use with samples that contain Tris buffer, heavy metals, or sulfides.



Fisher Scientific accumet AB200 pH/Conductivity/Temp Meter

An Accurate, Affordable, Multi-Parameter Instrument

- » New model measures pH & Conductivity/TDS/Resistivity/Salinity
- » Choose to display one parameter or pH and another parameter together
- » Graphic LCD with adjustable backlight
- » Date and time to meet GLP
- » Password protection
- » Optional calibration reminder & high/low alarm
- » 500 point memory – log manually or at intervals
- » Easily view calibration data

For best results, connect optional 13-620-BSP stirrer (see page 13) for efficient mixing and fast, accurate measurements!



Kit includes 13-620-100 cell.
For other conductivity cells, see page 34.

Specifications and Ordering Information

pH Mode	
Range	-2.000 to 20.000 pH
Resolution	0.1 / 0.01 / 0.001 pH
Accuracy	±0.002 pH + 1 LSD
Buffer Sets	USA: 2.000, 4.010, 6.997, 10.013, 12.000 NIST: 1.678, 4.010, 6.865, 9.184, 12.460 DIN: 1.090, 3.060, 4.650, 6.790, 9.230, 12.750 FSCI: 1.000, 3.000, 6.000, 8.000, 10.000, 13.000 CUSTOM: Any 2-5 values, ≥1.0 pH unit apart
Slope Display	Yes, up to 5 different slopes with offset
Temp Compensation	Automatic or manual (0 to 100 °C / 32 to 212 °F)
mV Mode	
Range	±2000.0 mV / Rel. mV
Resolution	0.1 mV
Accuracy	±0.2 mV or ±0.05 % whichever is greater
Offset Adjustment	Up to ±150 mV
Temperature Mode	
Resolution	0.1 °C / 0.1 °F
Accuracy	±0.3 °C / ±0.5 °F
Calibration	Offset in 0.1 ° increments; offset range: ±5 °C / 9 °F
Conductivity Mode	
Range	0.00 µS to 500.0 mS
Resolution	0.01 / 0.1 µS ; 0.001 / 0.01 / 0.1 mS
Accuracy	±1 % full scale
Coefficient (Per °C)	Linear & pure ; adjustable 0.000 to 10.000 %
Normalization	15.0 to 30.0 °C
TDS Mode	
Range	0.00 ppm to 500 ppt (@ TDS factor 1.00)
Resolution	0.01 / 0.1 ppm ; 0.001 / 0.01 / 0.1 ppt
Accuracy	±1 % full scale
TDS Factor	0.400 to 1.000
Salinity Mode	
Range	0 to 80.0 ppt
Resolution	0.01 / 0.1 ppm ; 0.001 / 0.01 / 0.1 ppt
Accuracy	±1 % full scale
Resistivity Mode	
Range	2.000 Ω to 20.0 MΩ
Resolution	0.01 / 0.1 Ω ; 0.001 / 0.1 kΩ ; 0.01 MΩ
Accuracy	±1 % full scale
Output	RS-232 (phono plug), mini-B USB, stirrer
Language Selection	English, Deutsch, 中文, Français, Italiano, & Español
Memory	500 data sets, viewable
Datalogging	Manual, timed (selectable every 3 to 3600 seconds) Printer or CSV format
Cal Due Alarm	Yes, user selectable from 8 hrs, 16 hrs, & 1-31 days
High / Low Alarms	Yes, user selectable, visual & audible

AB200 Meter Kit: Includes meter, TRIS compatible pH/ATC electrode (13-620-631), Conductivity/Temp probe (13-620-100), electrode arm, RS-232 & USB cables, 100/240 V power supply, and manual.

AB200 Meter Only: Includes meter, electrode arm, RS-232 & USB cables, 100/240 V power supply, and manual.

Description	Catalog No.
AB200 Kit	13-636-AB200
AB200 Meter Only	13-636-AB200A
Fisher Scientific accumet™ Benchtop Stirring Probe	13-620-BSP

The AB200 is a multi-parameter instrument capable of measuring pH, temperature, conductivity, total dissolved solids, resistivity, and salinity. Temperature is measured in two channels; a high accuracy thermistor built-in to

the conductivity probe, or separately along with the pH reading.

A generous graphic backlit display makes the AB200 meter easy to read, while plain language

prompts and error messages guide the user through measurements and calibrations. Change the display view with a press of a button.

Fisher Scientific accumet Benchtop Stirring Probe

Simple "Plug and Stir" Technology Available With New AB & XL Series

- » Provides mixing without magnetic stir bars or stir plates required!
- » Saves valuable bench space
- » Quiet operation
- » Adjustable speed control
- » No additional power required
- » Multi-parameter XL benchtop models can accept up to two probes for simultaneous stirring
- » Compatible with the following models:
 - XL150, XL200, XL250, XL500, XL600
 - AB150, AB250, AB200

Ordering Information

Description	Catalog No.
Fisher Scientific accumet Benchtop Stirring Probe	13-620-BSP
Replacement Paddle with Stirring Rod	13-620-RP
Electrode Holding Arm and Bracket for AB & XL Series Benchtop Meters	13-637-671



Slim design allows effortless mixing.





Fisher Scientific accumet AP110, AP115 And AP125 Meters

Accurate Measurements Anytime, Anywhere

- » Large LCD with backlighting and plain-language text – easy to view and use
- » Rugged waterproof IP67 housing
- » 0.001 pH resolution and date/time to meet GLP (AP115, AP125)
- » 200 data point internal memory
- » 3-year meter warranty
1-year electrode warranty



Specifications and Ordering Information

pH Mode	
Range	AP110 -2.00 to 20.00 AP115, AP125 -2.00 to 20.000
Resolution	AP110 0.1 / 0.01 AP115, AP125 0.1 / 0.01 / 0.001
Accuracy	AP110 ±0.01 AP115, AP125 ±0.002
Calibration	Up to 5 or 6 points from 3 sets: (USA: 2.00, 4.00, 7.00, 10.00 12.00; NIST: 1.68, 4.01, 6.86, 9.18, 12.45; EURO: 1.00, 3.00, 6.00, 8.00, 10.00, 13.00)
Ion Concentration Mode (AP125 Only)	
Range	0.001 to 99,999
Resolution	1, 2, or 3 significant digits
Accuracy	0.5 % FS (monovalent); 1 % FS (trivalent)
Calibration Points	From 2 to 5 points
Available Calibration Values	0.1, 0.5, 1, 2, 5, 10, 50, 100, 500, 1000
mV, Relative mV Mode	
Range	±2000
Resolution	0.1 / 1
Accuracy	±0.2 / 2
Calibration Points	1 (Rel mV)
Temperature Mode	
Range	-5 to 100 °C
Resolution	0.1 °C
Accuracy	±0.3
Calibration Points	1 (ATC) & 1 (MTC)
Available Calibration Values	1 point ±5 adjustment (ATC), -5 to 100 default adjustment (MTC)
Date/Time & GLP (AP115 & AP125 Only)	Yes

Meter Only: Includes meter, 9 V battery, and manual.

Meter Kit: Includes meter, pH/ATC electrode (13-620-AP50A), refill solution, hard carrying case, sample bottles, pH calibration packets, 9 V battery, and manual.

Description	Catalog No.
AP125 Meter Kit	13-636-AP125
AP125 Meter Only	13-636-AP125A
AP115 Meter Kit	13-636-AP115
AP115 Meter Only	13-636-AP115A
AP110 Meter Kit	13-636-AP110
AP110 Meter Only	13-636-AP110A

Replacement Parts And Accessories For AP100 Series Portable Meters

Description	Catalog No.
pH/ATC Electrode, Single-Junction, Epoxy Body, Refillable	13-620-AP50A
pH/ATC Electrode, Double-Junction, Epoxy Body, Refillable	13-620-AP61
pH/ATC Electrode, Double-Junction, Epoxy Body, Gel Filled	13-620-AP52
Optional 110/220 V Power Supply	13-636-100
Temperature Probe, Stainless Steel	13-620-AP53

Go Portable! Indoor or outdoor these new Fisher Scientific accumet meters are versatile and rugged enough for the most demanding applications. Soft touch keys with comfort grip make one-handed operation a breeze.

Meters are small enough for a shirt pocket, yet smart enough for any laboratory. The AP100 kits include an unbreakable pH/temperature electrode that is refillable for a fast response and the longest life.

The AP125 accepts any ion selective electrode (ISE) with a BNC connection for direct ion concentration measurement from 0.001 to 99,999 ppm.

Fisher Scientific accumet AP71 And AP72 Meters

Durable And Reliable Portable Meters



- » Measures pH, mV, and relative mV to 0.01 pH and 0.1 mV
- » Dual display shows pH or mV, and temperature
- » Push-button pH calibration with auto-buffer recognition
- » Manual or automatic temperature compensation
- » Store and recall 16 readings with corresponding temperature

AP72 Meter Also Features:

- » Real-time clock stamps stored data and calibration data with date and time
- » Expanded memory stores up to 50 readings
- » Selectable pH buffer sets: Standard US, NIST, and DIN buffers
- » Selectable automatic endpoint freezes reading when stable
- » °C/°F selectable

See pages 24 for a list of other pH electrodes.
See page 36 for calibration solutions.



- **Excellent ±0.01 pH accuracy**
- **Advanced setup mode allows for meter customization**
- **Accepts standard pH or ORP electrodes with a BNC connector**

The AP71 and AP72 meters feature advanced pH meter functions in a waterproof, dustproof

housing that floats. These meters read pH, mV or relative mV, and temperature in °C. The AP72 meter also reads temperature in °F. Both meters have excellent ±0.01 pH accuracy.

The ergonomic design of the AP71 and AP72 meter fits comfortably in your hand. Use the detachable electrode bracket to position the electrode for one-handed operation. Advanced setup mode lets

you customize meter parameters and check electrode slope and offset to ensure electrode accuracy.

The AP71 and AP72 pH meters are ideal for basic pH and mV readings in both harsh environmental conditions and general laboratory applications.

Specifications and Ordering Information

Memory	
AP71	Stores 16 data sets
AP72	Stores 50 data sets
pH Mode	
Range	-2.00 to 16.00
Resolution	0.01
Relative Accuracy	±0.01
Automatic Buffer Recognition	Yes
Calibration Points	
AP71	5 (1.68, 4.01, 7.01, 10.01, 12.45)
AP72	6 (US: 1.68, 4, 7, 10, 12.45; NIST: 1.68, 4.01, 6.86, 9.18, 12.45; or DIN: 1.09, 2.06, 4.65, 6.79, 9.23, 12.75)
mV Mode	
Range	±2000.0
Resolution	0.1 / 1
Accuracy	±0.2 / ±2
Temperature Mode	
Range	
AP71	0.0 to 100.0 °C
AP72	0.0 to 100.0 °C, 32.0 to 212.0 °F
Resolution	0.1 °C / 0.1 °F
Accuracy	±0.5 °C / ±0.5 °F
Inputs	BNC, ATC

Meter Only: Includes meter, AAA batteries, and manual.

Meter Kit: Includes meter, pH/ATC electrode (13-620-AP55), refill solution, hard carrying case, 60 ml sample bottles of storage solution, rinse water, pH 4 and pH 7 buffers, AAA batteries, and manual.

Description	Catalog No.
AP72 Meter Kit	13-636-AP72
AP72 Meter Only	13-636-AP72A
AP71 Meter Kit	13-636-AP71
AP71 Meter Only	13-636-AP71A

Replacement Parts And Accessories For AP 71 And AP 72 Series Portable Meters

Description	Catalog No.
Temperature Probe For AP 71 And AP 72, 316 Stainless Steel, 3-Ft Cable	13-620-20





Fisher Scientific accumet AP85 And AP75 Meters

Rugged Waterproof Portable Meters

- » Completely waterproof and dustproof
- » °C/°F selectable
- » ±1 full-scale accuracy for conductivity and TDS measurements
- » Switch from conductivity and TDS with the press of a button
- » Dual display shows conductivity or TDS plus temperature
- » Auto-ranging for conductivity across five ranges
- » Five point conductivity and TDS calibration

AP85 Meter Also Features:

- » pH measurement with five-point pH calibration and auto-buffer recognition

AP75 Meter Also Features:

- » Real-time clock stamps stored data and calibration data with date and time
- » Memory function stores and recalls up to 50 readings



AP85



AP85 meter kit

Specifications and Ordering Information

Memory (AP75 Only)	Stores 50 data sets
Conductivity Mode	
Cell Constants	
AP75	0.1, 1.0, 10
AP85	1.0
Range	
Conductivity	0.00 to 199.9 mS/cm
TDS	0.00 to 200 ppt
Accuracy	±1 % full scale
Calibration Points	Up to 5 points
Temperature Mode	
Range	0.0 to 100.0 °C, 32.0 to 212.0 °F
Resolution	0.1 °C / 0.1 °F
Accuracy	±0.5 °C / ±0.9 °F ±0.2 °C (±0.3 °F)
pH Mode (AP85 Only)	
Range	-2.00 to 16.00
Resolution	0.01
Relative Accuracy	±0.01
Calibration Points	5 (1.68, 4.01, 7.00, 10.00, 12.45)

AP85 Meter: Includes meter, pH/ATC electrode (13-620-AP55), conductivity/TDS/temp probe (13-620-AP54), AAA batteries, and manual.

AP85 Meter Kit: Includes meter, pH/ATC electrode (13-620-AP55), conductivity/TDS/temp probe (13-620-AP54), hard carrying case, pH buffer solutions, conductivity calibration solutions, AAA batteries, and manual.

AP75 Meter: Includes meter, conductivity/TDS/temp probe (13-620-AP54), batteries, and instruction manual.

AP75 Meter Kit: Includes meter, conductivity/TDS/temp probe (13-620-AP54), hard carrying case, conductivity calibration solutions, sample bottle, batteries, instruction manual, and hard plastic carrying case.

Description	Catalog No.
AP85 Meter Only	13-636-AP85A
AP85 Meter Kit	13-636-AP85
AP75 Meter Only	13-636-AP75A
AP75 Meter Kit	13-636-AP75

Replacement Parts And Accessories For AP85 And AP75 Meters

Description	Catalog No.
Replacement Conductivity/TDS/Temperature Probe	13-620-AP54
Replacement pH/ATC Electrode, 3-Ft Cable Electrode	13-620-AP55

The AP75 and AP85 meters measure both conductivity and TDS plus temperature with a housing that's perfect for rough conditions. These waterproof, dustproof meters float, and are very easy to clean.

Advanced setup mode lets you customize meter parameters. Auto-ranging for conductivity and TDS measurements offers fast response over the entire range. Adjust temperature coefficient from 0 to 10 % per °C or °F for accurate compensation in almost any solution.

The AP75 meter offers a real-time clock, expanded memory function, and selectable cell constant. The AP85 meter measures pH, conductivity and TDS plus temperature. AP85 also features five-point pH calibration with ATC and auto-buffer recognition.

Fisher Scientific accumet AP84 And AP74 Meters

Choose AP84 For pH/DO/Temp Or AP74 For DO/Temp Only



- » Durable Waterproof design with IP67 rating
- » No meter warm-up required
- » Adjustable backlit display
- » Dual display shows mg/L (ppm) or % saturation plus temperature
- » °C/°F selectable
- » Independent 100 % and zero adjustment calibrations
- » Key in salinity and pressure manually; the meter does the rest
- » 3-yr warranty

AP84 Meter Also Features:

- » pH measurement with auto-buffer recognition
- » Meter displays probe diagnostics, view slope, offset, and mV values



AP84

AP74 Meter Also Features:

- » Real-time clock stamps stored data and calibration data with date and time
- » Memory function stores and recalls up to 50 readings



AP74

The AP74 and AP84 meters read dissolved oxygen in mg/L, ppm, or % saturation. They use barometric pressure and salinity offset to ensure high accuracy with independent 100 %, zero and offset adjustment capabilities. The adjustable backlit display is ideal for taking readings in low light. Probe includes 10-ft standard cable length.

These meters also feature a galvanic probe design. This design allows the user to take dissolved oxygen measurements immediately, without the warm up time of other probe designs! Advanced features include display of probe diagnostics. The waterproof, dustproof housing floats, and is very easy to clean.

The AP84 meter also takes pH readings with ± 0.01 pH accuracy.

Specifications and Ordering Information

Memory (AP74 Only)	Stores 50 data sets
Dissolved Oxygen Mode	
Range	
mg/L (ppm)	0.00 to 20.00
% Saturation	0.0 to 200.0
Resolution	
mg/L (ppm)	0.01
% Saturation	0.1
Accuracy	± 1.5 % full-scale
Calibration Points	2 point (0 %, 100 %), 1 point (mg/L)
Temperature Mode	
Range	0.0 to 50.0 °C (32.0 to 122.0 °F)
Resolution	0.1 °C (0.1 °F)
Accuracy	± 0.3 °C (± 0.5 °F)
pH Mode (AP84 Only)	
Range	-2.00 to 16.00
Resolution	0.01
Relative Accuracy	± 0.01
Calibration Points	Up to 5 (1.68, 4.01, 7.00, 10.00, 12.45)

Meter: Includes meter, dissolved oxygen/temp probe (15-500-034), electrolyte, (2) membrane cap assemblies (15-500-037), polishing disk, AAA batteries, and manual.

Meter Kit: Includes meter, dissolved oxygen/temp probe (15-500-034), pH electrode (AP84 only) (13-620-AP56), electrolyte, (2) membrane cap assemblies (15-500-037), hard carrying case, polishing disk, AAA batteries, and manual.

Description	Catalog No.
AP84 Meter Kit	13-636-AP84
AP84 Meter	13-636-AP84A
AP74 Meter Kit	13-636-AP74
AP74 Meter	13-636-AP74A

Replacement Parts And Accessories For AP74 And AP84 Meters

Description	Catalog No.
Replacement Dissolved Oxygen/Temperature Probe With 10-Ft Submersible Cable	15-500-034
Electrolyte Solution For Dissolved Oxygen Probes, 500 mL	15-500-036
Replacement pH Electrode For AP84, ABS Body, 10-Ft Cable, 25 mm Dia.	13-620-AP56



AP84 meter kit

Fisher Scientific accumet Electrodes

pH Electrodes For Virtually Any Sample Type

When selecting a pH system, choose your meter based on what features you need, i.e. resolution, output, memory, etc. Choose the corresponding pH electrode for your meter based primarily on your sample type and conditions, i.e. wastewater with sulfides, room temperature, student use, 5 days/week, etc.

Take a glance in the Fisher Scientific catalog and you will find 100's of pH electrodes to choose from. While many electrodes might be work adequately for a particular application, not all will perform equally or last as long as others. Usually in situations in which a pH electrode "didn't last long", the electrode is not matched well for the application resulting in poor performance, and ultimately failure. Understanding the different electrode options that are available and knowing how to use them to your advantages is a critical step to getting the most out of your pH measurement system.


The following guide is designed to help you understand the ABC's of electrode selection. If you are still undecided or have questions regarding any product, your Fisher Scientific representative and our electrochemistry experts (888-358-4706 or accumet@fishersci.com) are there to help guide you!

» The ABC's Of Electrode Selection

Combination Or Half-Cell

There are two components within a pH electrode system. A pH indicating (or sensing) electrode develops a potential dependent on the pH, and the reference electrode which provides a constant potential to completes the electrical circuit. Combination electrodes have both the indicating and reference electrodes "combined" into one electrode. Alternatively, separate half-cell electrodes can be used. Since reference electrodes often outlast sensing electrodes, replacing indicating electrodes can mean lower replacement costs than replacing an equivalent combination electrode.

In practical terms, nearly all electrodes used today are combination electrodes. There are many reasons including; many pH meters require an adapter to accept half-cell electrodes, handling multiple electrodes is un-desirable (or impossible with small samples!), half cells don't have built in ATC and therefore require a third electrode for temperature compensation, the complexity of diagnosing electrode problems and most of all, the reduced cost and performance of today's combination electrode designs.

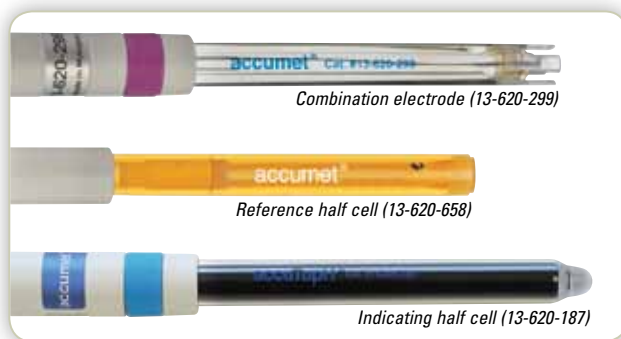

Verdict: Use a combination electrode unless the method you must follow calls for half-cells. You'll have many more choices available to you. Combination cells may or may not have a temperature sensor built-in.

Glass Or Plastic Body

It probably goes without saying, but if an electrode literally breaks into pieces, it is useless and can not be repaired. Combination glass and combination plastic electrodes use an indicating electrode with a glass sensing bulb on the end. This is important for several reasons. First, plastic electrodes are not immune from breakage. Second, if an electrode breaks it will likely be at the tip, not the body itself. A plastic electrode with little to no bulb protection defeats the purpose of a plastic electrode in the first place.

To decide on which to use, let us look at the advantages and disadvantages of each, starting with the glass body electrode. Glass electrodes are easier to clean and maintain since they can tolerate just about any solvent and inorganic material (with the exception of HF!) and can handle higher temperatures quite nicely – typically to 100 °C. The fact that glass electrodes also have a glass sensing bulb is also an advantage. Since the seal that combines the bulb to the body is similar material, it is one less thing that can go wrong during measurement and doesn't become the source of junction potential as it does in plastic electrodes. This is especially important consideration for applications that have repeated and extreme heating and cooling – the expansion and contraction that occurs is handled much better by glass electrodes. The downside of glass electrodes is fairly easy – they are generally more expensive than plastic, and they have a greater potential for breakage.

Plastic electrodes are less expensive than glass equivalents and can usually take much abuse in the lab and in the field. Most electrodes with built-in temperature compensation elements are plastic due to the complexity in manufacturing them. As a result, they are most popular with field and portable meters, but can also be used in laboratory environments (such as 13-620-631). To protect the glass sensing bulb, many plastic electrodes use an integral housing that limit the bulb exposure, but often can be difficult to clean.

Verdict: Glass electrodes are definitely worth the upgrade if you have significant temperature fluctuations. If bulb breakage is a concern, consider Fisher Scientific accumet accuTupH electrodes with thick glass bulbs! If you want ATC built-in to your electrode, expect to settle for plastic.

Refillable Or Non-Refillable (Gel)

All pH electrodes use/leak solution. Refillable electrodes do so more quickly, and can be replenished when they require more filling solution. Gel filled electrodes do so very slowly and when they run out or the gel is no longer flowing, can not be replenished and must be replaced.

Refillable electrodes are generally more expensive than gel-filled equivalent electrodes but respond much faster. They also last longer, because the filling solution can be replaced indefinitely; however the periodic addition of filling solution that is required also happens to be the main disadvantage. Another downside is that when the filling hole is left open for an extended period, dried salt may be left behind which often involves cleaning. The act of refilling and opening and closing the fill hole with Fisher Scientific accumet electrodes is extremely easy due to the patented filling mechanism. It takes just seconds to open the hole and a few seconds more to fill the probe.

Gel-filled electrodes are less expensive, require less maintenance, and are usually plastic. High quality gel formulations have also extended the once limited shelf-life in recent years.

Verdict: Refillable electrodes are usually worth the extra maintenance – especially if it's a Fisher Scientific accumet electrode.

Fisher Scientific accumet Electrodes

pH, pH/ATC, ORP, and Ion Selective Electrodes


Single Or Double-Junction (Tris Compatible)



The single-junction electrode on the left has a black, clogged junction and is no longer responsive.

This decision is extremely important and should not be overlooked. If you will be measuring samples that have sulfides, proteins, heavy metals, TRIS, or anything that might react with silver, or if you will be testing samples that are unknown, use a double-junction electrode. Calomel electrodes would also be suitable but have fallen out of favor due to mercury content and regulations that ban shipments of them in specific states in the US. Single-junction electrodes are less expensive, but offer no other advantages. If you use a single-junction electrode in a solution with TRIS, it's just a matter of time before it fails.

Verdict: If you will only measure drinking water, you can save money by using a single-junction pH electrode. If you have TRIS, sulfides, proteins, heavy metals or are measuring samples that are unknown, look for a Fisher Scientific accumet electrode with a purple ring – indicating that is it compatible.

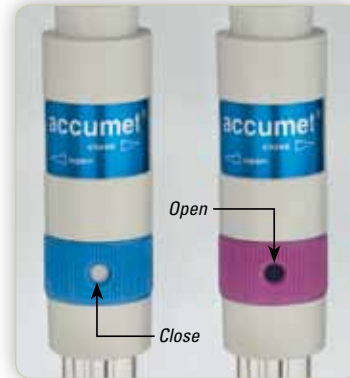

Refillable electrodes use our patented twist open and close mechanism.

Color coded electrode bands simplify electrode selection:

- ◆ Purple = TRIS Compatible
- ◆ Blue = General Purpose



Over 30 years of experience in the design, development, and manufacture of electrodes go into each Fisher Scientific accumet electrode.

We offer electrodes that provide fast, accurate measurements in hundreds of different applications – including yours!

A complete line for every application: made with care and precision. All Fisher Scientific accumet electrodes feature continuous electrical shielding and insulation of the internal elements, cable and connectors for extremely stable, reproducible readings with a minimum of electrical noise. Each electrode is individually tested, serialized to meet GLP requirements, and backed by a knowledgeable support staff (888-358-4706 or accumet@fishersci.com) and 1 year warranty.





Fisher Scientific accuPest pH Electrodes

High Performance Models For Critical Research

State-of-the-art design for fast, accurate measurements despite sample temperature differences – plus extra durability. Feature innovative reference system that controls chemical equilibria, prevents precipitation of solution components at reference element from 0 to 100 °C; plus internal electrolyte with minimal temperature coefficient. Result: highly predictable, super reliable electrodes that respond quickly at any temperature. Cycle between 25 and 80 °C samples, reach reproducible pH in 30 seconds (vs. 1 to 3 minutes

for other electrodes). Drift and accuracy problems are virtually eliminated.

Read sample pH in <20 seconds, correct to ±0.02 pH; pH value stays constant at any temperature. Best of all, these electrodes read pH consistently at elevated temperatures – and without premature loss in performance.

Choice of standard-size glass body, epoxy body with flushable junction, and glass body with flushable junction.

accuPest R pH Electrodes

Catalog No.	13-620-195	13-620-196	13-620-197
Special	accuPest R	accuPest R accuFlow	accuPest R accuFlow
Parameter	pH	pH	pH
Combination Or Half Cell	Combination	Combination	Combination
ATC Connection	n/a	n/a	n/a
Refillable Or Gel (Sealed)	Refillable	Refillable	Refillable
Refill Solution	SP138-500	SP138-500	SP138-500
Junction Type	Double	Double	Double
Glass Or Plastic Body	Glass	Plastic	Glass
Max Temp	100 °C	80 °C	100 °C
Length x Diameter (mm)	102 x 12	102 x 12	102 x 12
Note:	High performance, ideal for samples temp variation	High performance, flushable junction for tough samples	High performance, flushable junction for tough samples

Fast and accurate for samples at widely varying temperatures. Patented design: dual ceramic junctions, sealed reference, and special internal electrolyte to eliminate slow response when measuring samples at different temperatures in quick succession. Accurate to ±0.01 pH at 25 °C and ±0.05 pH from -5 to 100 °C. Response times of 20 seconds or less. Negligible drift.

Isolated reference and outer KCl fill solution prevent clogging from silver-compound precipitates. Unique pH bulb is filled with special crystals to speed thermal equilibrium. Choice of four styles: standard-size glass body, MicroProbe™ extra-long glass body, extra-long epoxy body, and pH/ATC epoxy body.

accuPest pH Electrodes

Catalog No.	13-620-296	13-620-297	13-620-298	13-620-113	13-620-114
Special	accuPest	accuPest long & narrow	accuPest long	accuPest	accuPest
Parameter	pH	pH	pH	pH/ATC	pH/ATC
Combination Or Half Cell	Combination	Combination	Combination	Combination	Combination
ATC Connection	n/a	n/a	n/a	13-620-16	13-620-19
Refillable Or Gel (Sealed)	Refillable	Refillable	Refillable	Refillable	Refillable
Refill Solution	SP138-500	SP138-500	SP138-500	SP138-500	SP138-500
Junction Type	accuPest (Tris compatible)	accuPest (Tris compatible)	accuPest (Tris compatible)	accuPest (Tris compatible)	accuPest (Tris compatible)
Glass Or Plastic Body	Glass	Glass	Plastic	Plastic	Plastic
Max Temp	100 °C	80 °C	80 °C	80 °C	80 °C
Length x Diameter (mm)	102 x 10	165(L) x 75 x 5	140 x 10	143 x 10	143 x 10
Note:	High performance, ideal for samples temp variation	High performance	High performance	See page 34 for list of discontinued meters using 13-620-16 ATC	ATC fits XL, AB, and AR meters

Fisher Scientific accumet pH Electrodes

Top Selling Rugged Glass And Capillary Junction Electrodes



Five times thicker than conventional glass pH electrodes. For applications where glass bulbs break frequently and epoxy body electrodes aren't practical. Up to 40 times tougher than conventional glass pH electrodes, without sacrificing response times.

13-620-185 also utilizes accupHast temp reference for top performance from a rugged glass electrode.


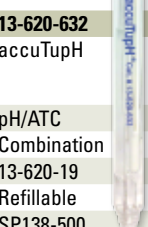





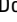





8-mil standard electrode



40-mil accuTupH rugged bulb

accuTupH+ & accuTupH pH Electrodes

						
Catalog No.	13-620-185	13-620-632	13-620-183A	13-620-182A	13-620-181	13-620-187
Special	accuTupH+, accupHast	accuTupH	accuTupH	accuTupH, US Standard Connector	accuTupH	accuTupH
Parameter	pH	pH/ATC	pH	pH/ATC	pH	pH indicating
Combination Or Half Cell	Combination	Combination	Combination	Combination	Combination	Half Cell, BNC
ATC Connection	n/a	13-620-19	n/a	n/a	n/a	n/a
Refillable Or Gel (Sealed)	Refillable	Refillable	Refillable	Refillable	Refillable	n/a
Refill Solution	SP138-500	SP138-500	SP138-500	SP138-500	SP135-500	n/a
Junction Type	accupHast (Tris compatible) 	Double 	Double 	Double 	Single 	n/a
Glass Or Plastic Body	Glass	Glass	Glass	Glass	Glass	Glass
Max Temp	100 °C	100 °C	100 °C	100 °C	100 °C	100 °C
Length x Diameter (mm)	102 x 10	102 x 12	102 x 12	102 x 10	102 x 10	106 x 12
Note:	Rugged & fast temp response, high performance, includes bulb protector	Rugged, Tris compatible, includes bulb protector. ATC for AB, AR, and XL meters.	Rugged, Tris compatible, includes bulb protector. Included with AB150B and XL kits	Rugged, Tris compatible, includes bulb protector. US Std Connector (not BNC)	Rugged, general purpose. includes bulb protector	Rugged, use with reference half cell








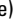
Single-pore capillary junction provides a flow channel about 200 times larger than typical reference junctions. Combined with a specially formulated

flowing gel reference electrolyte (13-636-430), provides a fast, virtually clog-free reference. The result is a faster, more stable pH measurement.

accuCap Capillary Junction Electrodes

Electrolyte & syringe (13-636-430)





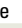
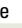

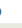
				
Catalog No.	13-620-130	13-620-131	13-620-132	13-620-133
Special	accuCap	accuCap	accuCap	accuCap, spear tip
Parameter	pH	pH	pH	pH
Combination Or Half Cell	Combination	Combination	Combination	Combination
ATC Connection	n/a	n/a	n/a	n/a
Refillable Or Gel (Sealed)	Refillable	Gel	Gel	Gel
Refill Solution	13-636-430	n/a	n/a	n/a
Junction Type	Capillary open pore (Tris compatible) 	Capillary open pore (Tris compatible) 	Capillary open pore (Tris compatible) 	Capillary open pore (Tris compatible) 
Glass Or Plastic Body	Glass	Glass	Plastic	Glass
Max Temp	80 °C	80 °C	60 °C	50 °C
Length x Diameter (mm)	160 x 12	130 x 12	120 x 12	80(L) x 25 x 6
Note:	Research quality, included with XL series pH kits	Non-refillable glass electrodes like this are hard to find	General purpose	Spear tip and 6 mm diameter useful for semi solids & small samples

Fisher Scientific accuTmet pH Electrodes


pH Electrodes To Match Your Application Type



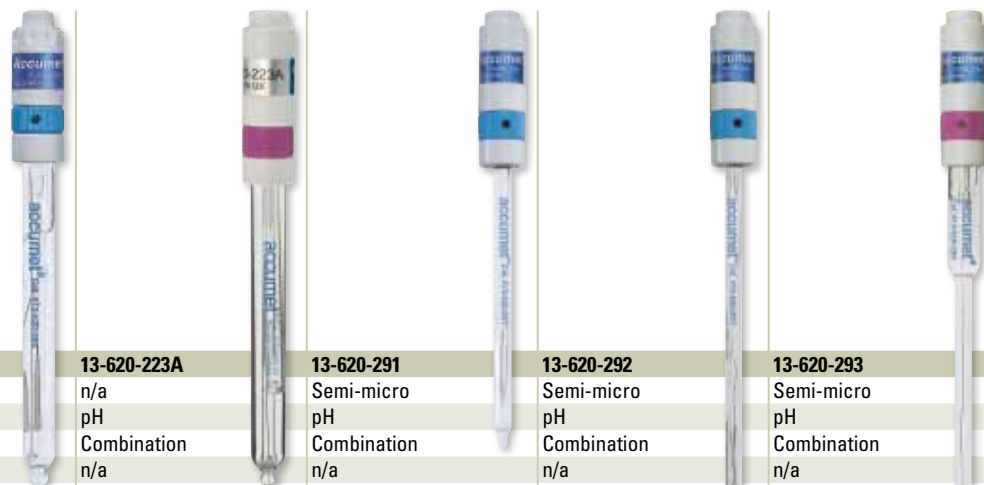


Refillable Glass pH Electrodes

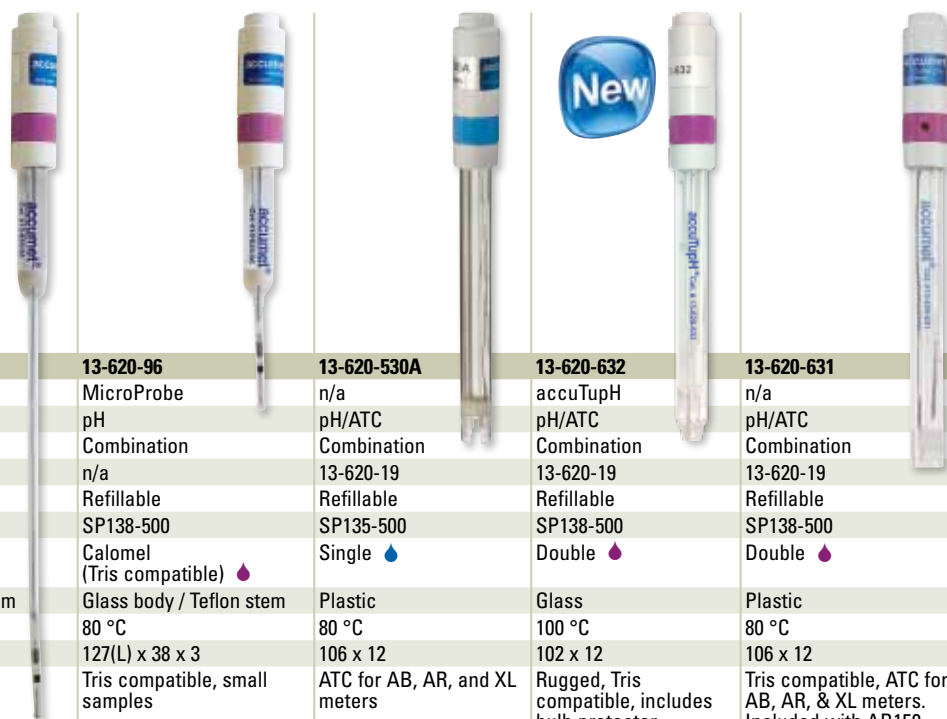


Catalog No.	13-620-285	13-620-223A	13-620-291	13-620-292	13-620-293
Special	n/a	n/a	Semi-micro	Semi-micro	Semi-micro
Parameter	pH	pH	pH	pH	pH
Combination Or Half Cell	Combination	Combination	Combination	Combination	Combination
ATC Type	n/a	n/a	n/a	n/a	n/a
Refillable Or Gel (Sealed)	Refillable	Refillable	Refillable	Refillable	Refillable
Refill Solution	SP135-500	SP138-500	SP135-500	SP135-500	SP138-500
Junction Type	Single 	Double 	Single 	Single 	Calomel 
Glass Or Plastic Body	Glass	Glass	Glass	Glass	Glass
Max Temp	100 °C	100 °C	100 °C	100 °C	80 °C
Length x Diameter (mm)	102 x 12	102 x 12	100 x 6	150 x 6	160(L) x 120 x 6
Note:	General purpose. Included with AB15+ and AB15 kits. Includes bulb protector	Tris compatible, includes bulb protector	6 mm diameter for small samples, test tubes	Same as 13-620-291 but longer	Tris compatible

Micro pH And pH/ATC Refillable Electrodes




Catalog No.	13-620-95	13-620-96	13-620-530A	13-620-632	13-620-631
Special	MicroProbe	MicroProbe	n/a	accuTupH	n/a
Parameter	pH	pH	pH/ATC	pH/ATC	pH/ATC
Combination Or Half Cell	Combination	Combination	Combination	Combination	Combination
ATC Type	n/a	n/a	13-620-19	13-620-19	13-620-19
Refillable Or Gel (Sealed)	Refillable	Refillable	Refillable	Refillable	Refillable
Refill Solution	SP138-500	SP138-500	SP135-500	SP138-500	SP138-500
Junction Type	Calomel (Tris compatible) 	Calomel (Tris compatible) 	Single 	Double 	Double 
Glass Or Plastic Body	Glass body / Teflon stem	Glass body / Teflon stem	Plastic	Glass	Plastic
Max Temp	80 °C	80 °C	80 °C	100 °C	80 °C
Length x Diameter (mm)	254(L) x 150 x 3	127(L) x 38 x 3	106 x 12	102 x 12	106 x 12
Note:	Tris compatible, small samples, long test tubes / NMR	Tris compatible, small samples	ATC for AB, AR, and XL meters	Rugged, Tris compatible, includes bulb protector. ATC for AB, AR, and XL meters.	Tris compatible, ATC for AB, AR, & XL meters. Included with AB150, AB200, & AB250 kits.

Refillable Plastic pH Electrodes

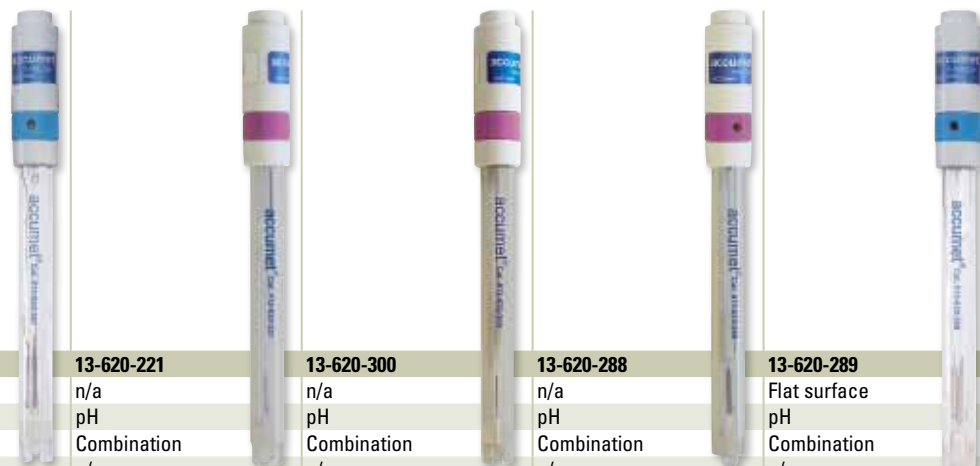

Catalog No.	13-620-287A	13-620-221	13-620-300	13-620-288	13-620-289
Special	n/a	n/a	n/a	n/a	Flat surface
Parameter	pH	pH	pH	pH	pH
Combination Or Half Cell	Combination	Combination	Combination	Combination	Combination
ATC Type	n/a	n/a	n/a	n/a	n/a
Refillable Or Gel (Sealed)	Refillable	Refillable	Refillable	Refillable	Refillable
Refill Solution	SP135-500	SP138-500	SP138-500	SP138-500	SP135-500
Junction Type	Single ◆	Double ◆	Calomel ◆	Calomel ◆	Single ◆
Glass Or Plastic Body	Plastic	Plastic	Plastic	Plastic	Plastic
Max Temp	80 °C	80 °C	80 °C	80 °C	80 °C
Length x Diameter (mm)	106 x 12	102 x 10	106 x 12	106 x 12	114 x 13
Note:	Same as 13-620-530A without ATC	Tris compatible, includes bulb protector	Tris compatible, with integral bulb guard	Tris compatible	Flat surface for agar, cheese, food, and more

Gel-Filled Plastic pH Electrodes

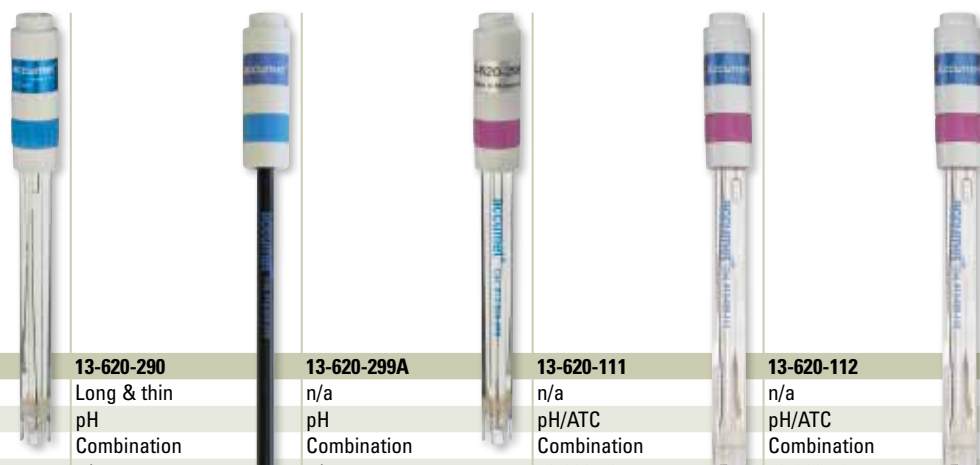

Catalog No.	13-620-108A	13-620-290	13-620-299A	13-620-111	13-620-112
Special	n/a	Long & thin	n/a	n/a	n/a
Parameter	pH	pH	pH	pH/ATC	pH/ATC
Combination Or Half Cell	Combination	Combination	Combination	Combination	Combination
ATC Type	n/a	n/a	n/a	13-620-19	13-620-16
Refillable Or Gel (Sealed)	Gel	Gel	Gel	Gel	Gel
Refill Solution	n/a	n/a	n/a	n/a	n/a
Junction Type	Single ◆	Single ◆	Double ◆	Double ◆	Double ◆
Glass Or Plastic Body	Plastic	Plastic	Plastic	Plastic	Plastic
Max Temp	80 °C	80 °C	80 °C	80 °C	80 °C
Length x Diameter (mm)	106 x 12	178 x 6	106 x 12	106 x 12	106 x 12
Note:	Economical, general purpose	Tall flasks, bottles	Economical, Tris compatible	ATC for AB, AR, and XL meters	See page 34 for list of discontinued meters using 13-620-16 ATC

Fisher Scientific accumet pH Electrodes





New pH/ATC Electrodes For Non-accumet Meters


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
- Combination pH mercury-free electrodes with built-in temperature compensation
- Fast, accurate response from 5 to 80 °C
- Double-junction pH/ATC electrodes compatible with Tris, proteins and sulfides

- Epoxy body is impact resistant and ideal for rough handling

All electrodes have a BNC connector and ATC connector; ATC will differ with meter type. 3-ft cable and electrode storage bottle are included, refillable models also include a 30 mL bottle of filling solution.

Universal pH/ATC Electrodes

Catalog No.	13-620-31C	13-621-701	13-621-702	13-621-703
Parameter	pH/ATC	pH/ATC	pH/ATC	pH/ATC
Combination Or Half Cell	Combination	Combination	Combination	Combination
ATC Connection	RCA (Cinch) plug	RCA (Cinch) plug	3.5 audio plug	Banana plug
Refillable Or Gel (Sealed)	Refillable	Gel	Gel	Gel
Refill Solution	SP135-500	n/a	n/a	n/a
Junction Type	Single 	Double 	Double 	Double 
Glass Or Plastic Body	Plastic	Plastic	Plastic	Plastic
Max Temp	80 °C	80 °C	80 °C	80 °C
Length x Diameter (mm)	106 x 12	106 x 12	106 x 12	106 x 12
Note:	ATC for Mettler™, Pinnacle™ and Corning™ meters	ATC for Mettler™, Pinnacle™ and Corning™ meters	ATC for Beckman™ meters	ATC for WTW™ and Pinnacle™ (part numbers ending with "P")


Catalog No.	13-620-111T	13-620-111	13-620-631
Parameter	pH/ATC	pH/ATC	pH/ATC
Combination Or Half Cell	Combination	Combination	Combination
ATC Connection	Mini DIN	Mini phone	Mini phone
Refillable Or Gel (Sealed)	Gel	Gel	Refillable
Refill Solution	n/a	n/a	SP138-500
Junction Type	Double 	Double 	Double 
Glass Or Plastic Body	Plastic	Plastic	Plastic
Max Temp	80 °C	80 °C	80 °C
Length x Diameter (mm)	106 x 12	106 x 12	106 x 12
Note:	ATC for Thermo Scientific Star meters	ATC for Denver / Sartorius meters	ATC for Denver / Sartorius meters





Fisher Scientific accumet pH And Half Cell Electrodes

pH/ATC Electrodes For Portable Fisher Scientific accumet Meters

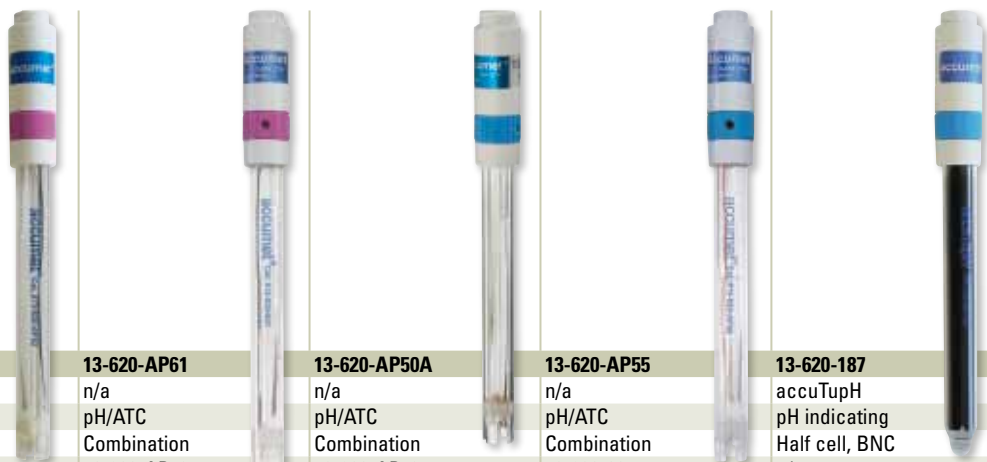




pH/ATC Electrodes For Field Use, accuTupH Half Cell



Catalog No.	13-620-AP52	13-620-AP61	13-620-AP50A	13-620-AP55	13-620-187
Special	n/a	n/a	n/a	n/a	accuTupH
Parameter	pH/ATC	pH/ATC	pH/ATC	pH/ATC	pH indicating
Combination Or Half Cell	Combination	Combination	Combination	Combination	Half cell, BNC
ATC Type	13-620-16	13-620-AP53	13-620-AP53	13-620-20	n/a
Refillable Or Gel (Sealed)	Gel	Refillable	Refillable	Refillable	n/a
Refill Solution	n/a	SP138-500	SP135-500	SP135-500	n/a
Junction Type	Double 	Double 	Single 	Single 	n/a
Glass Or Plastic Body	Plastic	Plastic	Plastic	Plastic	Glass
Max Temp	80 °C	80 °C	80 °C	80 °C	100 °C
Length x Diameter (mm)	102 x 12	102 x 12	102 x 12	102 x 12	106 x 12
Note:	ATC for AP60 & AP100 series meters	ATC for AP60 & AP100 series meters. Replaces 13-620-AP51	ATC for AP60 & AP100 series meters	ATC for AP70 & AP80 series meters	Rugged, use with reference half cell

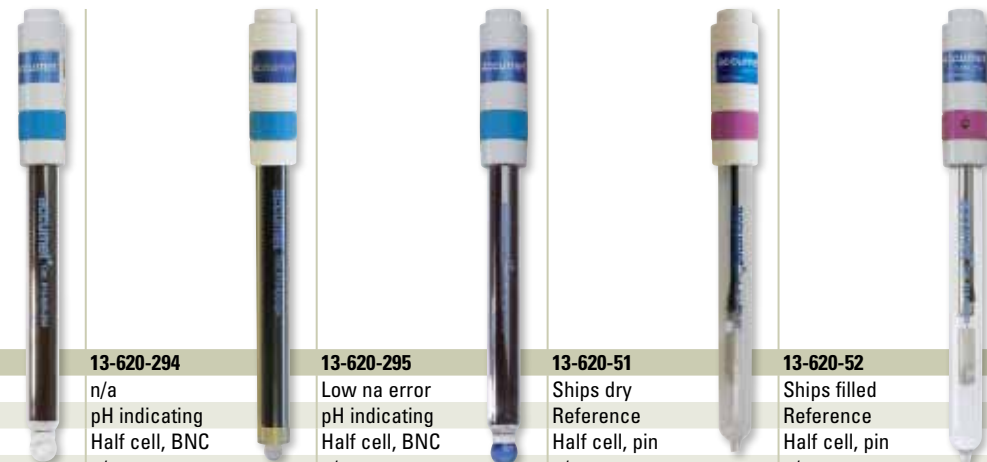Half-Cell Electrodes




Catalog No.	13-620-284	13-620-294	13-620-295	13-620-51	13-620-52
Special	n/a	n/a	Low na error	Ships dry	Ships filled
Parameter	pH indicating	pH indicating	pH indicating	Reference	Reference
Combination Or Half Cell	Half cell, BNC	Half cell, BNC	Half cell, BNC	Half cell, pin	Half cell, pin
ATC Type	n/a	n/a	n/a	n/a	n/a
Refillable Or Gel (Sealed)	n/a	n/a	n/a	Refillable	Refillable
Refill Solution	n/a	n/a	n/a	SP138-500	SP138-500
Junction Type	n/a	n/a	n/a	Calomel 	Calomel 
Glass Or Plastic Body	Glass	Plastic	Plastic	Glass	Glass
Max Temp	100 °C	80 °C	80 °C	80 °C	80 °C
Length x Diameter (mm)	106 x 12	106 x 12	106 x 12	106 x 12	106 x 12
Note:	Use with reference half cell	Use with reference half cell	Ideal for samples >pH 11, use with reference half cell	Pin connector, use with indicating half cell	Pin connector, use with indicating half cell



Fisher Scientific accumet Electrodes


Reference Electrodes Use Common Pin Connector Type

Half Cell Reference Electrodes

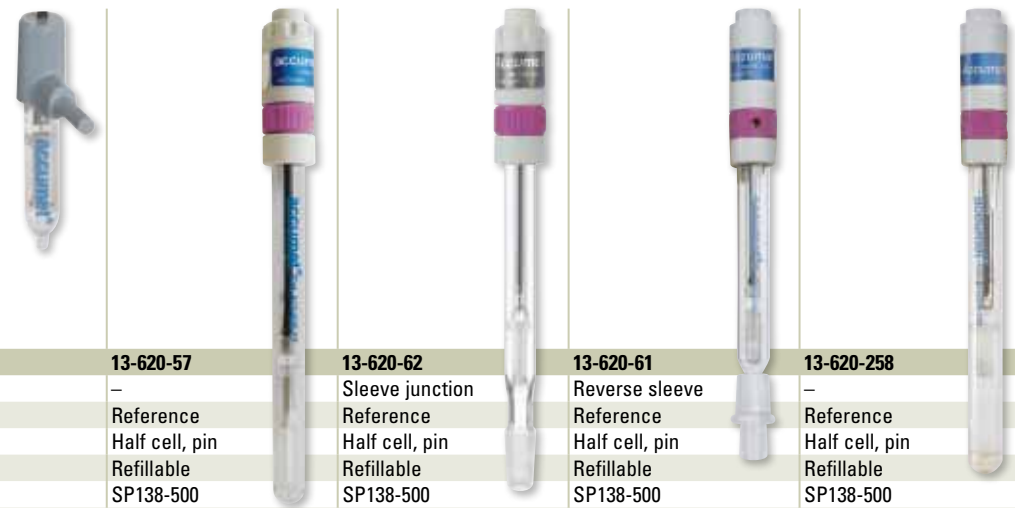

Catalog No.	13-620-79	13-620-57	13-620-62	13-620-61	13-620-258
Special	Minature	–	Sleeve junction	Reverse sleeve	–
Parameter	Reference	Reference	Reference	Reference	Reference
Combination Or Half Cell	Half cell, pin	Half cell, pin	Half cell, pin	Half cell, pin	Half cell, pin
Refillable Or Gel (Sealed)	Refillable	Refillable	Refillable	Refillable	Refillable
Refill Solution	SP138-500	SP138-500	SP138-500	SP138-500	SP138-500
Junction Type	Calomel	Calomel	Calomel	Calomel	Calomel
Glass Or Plastic Body	Glass	Glass	Glass	Glass	Plastic
Max Temp	80 °C	80 °C	80 °C	80 °C	80 °C
Length x Diameter (mm)	41 x 12	106 x 12	106 x 12, 16 (with sleeve)	106 x 12, 16 (with sleeve)	106 x 12
Note:	Pin connector, use with indicating half cell	Pin connector, non-aqueous samples, use with indicating half cell	Pin connector, for viscous samples, use with indicating half cell	Pin connector, for viscous samples, use with indicating half cell	Pin connector, use with indicating half cell


Half Cell Reference Electrodes

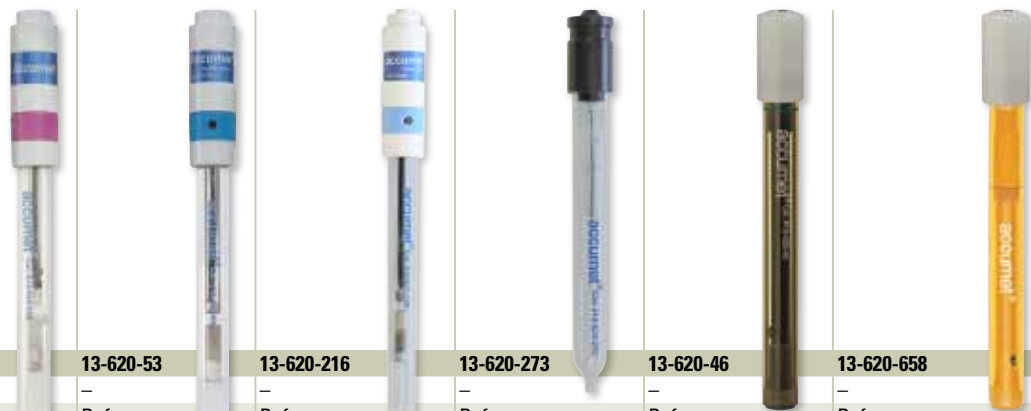

Catalog No.	13-620-259	13-620-53	13-620-216	13-620-273	13-620-46	13-620-658
Special	–	–	–	–	–	–
Parameter	Reference	Reference	Reference	Reference	Reference	Reference
Combination Or Half Cell	Half cell, pin	Half cell, pin	Half cell, pin	Half cell, pin	Half cell, pin	Half cell, pin
Refillable Or Gel (Sealed)	Gel	Refillable	Refillable	Refillable	Refillable	Refillable
Refill Solution	n/a	SP135-500	1M sodium sulfate	SP138-500	SP135-500	Inner 13-620-433 Outer 13-620-434
Junction Type	Calomel	Single	Single	Double	Single	Double
Glass Or Plastic Body	Plastic	Glass	Glass	Glass	Plastic	Plastic
Max Temp	80 °C	100 °C	100 °C	100 °C	100 °C	100 °C
Length x Diameter (mm)	106 x 12	106 x 12	106 x 12	106 x 12	108 x 13	108 x 13
Note:	Pin connector, use with indicating half cell	Pin connector, use with indicating half cell	Pin connector, use with indicating half cell (13-620-122 recommended) Ag and Cl titrations.	Pin connector, use with indicating half cell	Pin connector, use with indicating half cell	Pin connector, bromide, chloride, copper, iodide, lead, nitrate, silver/sulfide, redox, and pH applications requiring sample-compatible electrolyte, use with indicating half cell

Fisher Scientific accumet Half Cell And ORP Electrodes


Electrodes For Titrations, Redox, And Specialty Applications

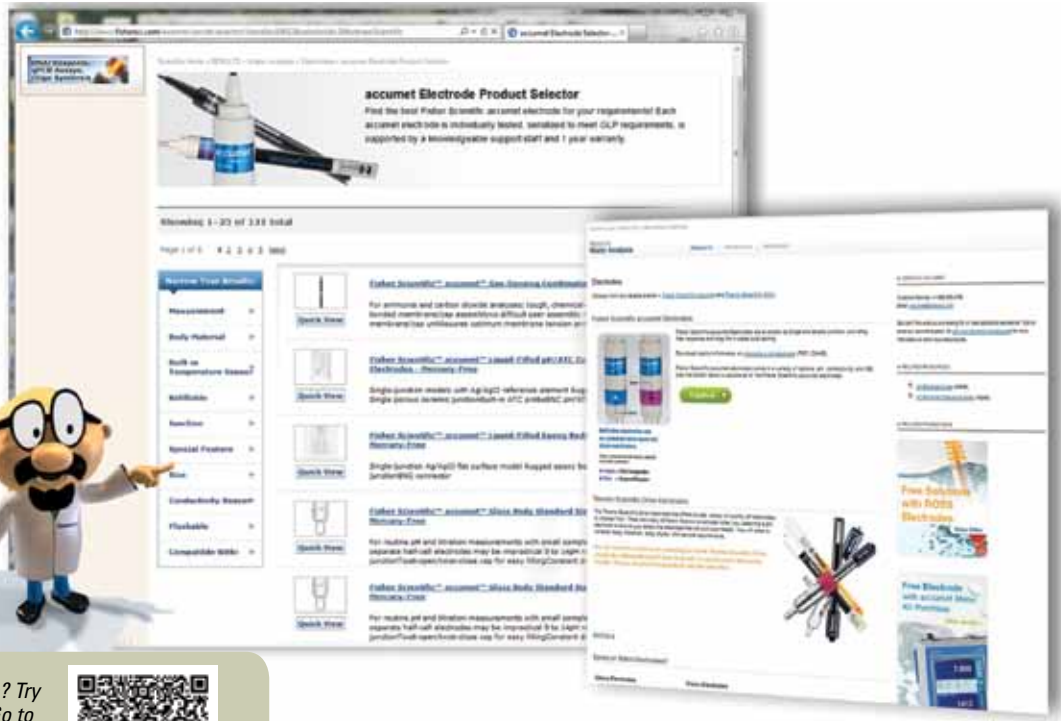

Trouble finding an electrode? Try our product selector tool. Go to fishersci.com/wateranalysis then click products, then electrodes. Or, scan this:



Reference, Metallic, And ORP Electrodes

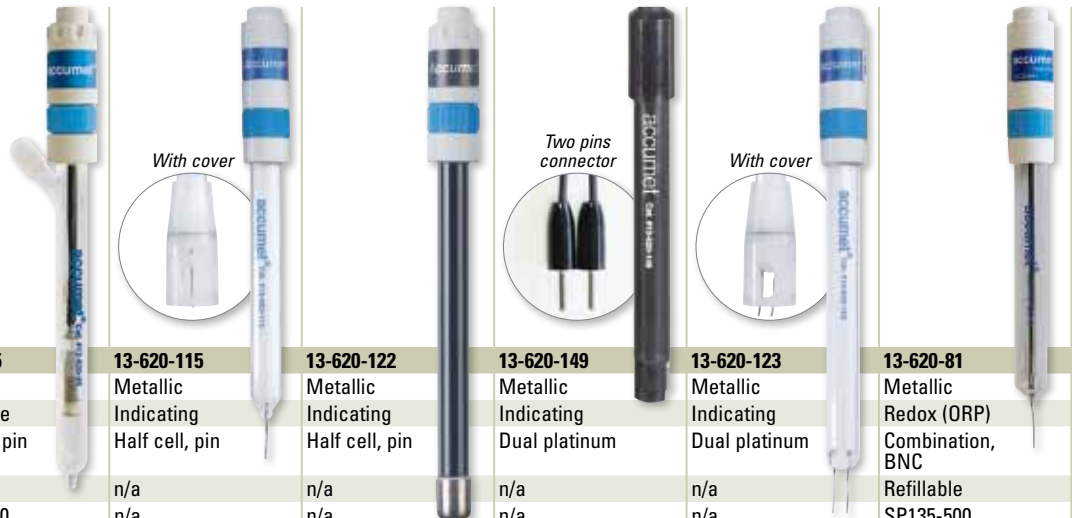

Catalog No.	13-620-45	13-620-115	13-620-122	13-620-149	13-620-123	13-620-81
Special	Side-arm	Metallic	Metallic	Metallic	Metallic	Metallic
Parameter	Reference	Indicating	Indicating	Indicating	Indicating	Redox (ORP)
Combination Or Half Cell	Half cell, pin	Half cell, pin	Half cell, pin	Dual platinum	Dual platinum	Combination, BNC
Refillable Or Gel (Sealed)	Refillable	n/a	n/a	n/a	n/a	Refillable
Refill Solution	SP135-500	n/a	n/a	n/a	n/a	SP135-500
Junction Type	Single	Redox	Redox	Redox	Redox	Redox
Glass Or Plastic Body	Glass	Glass	Glass	Plastic	Glass	Glass
Max Temp	100 °C	100 °C	100 °C	80 °C	100 °C	100 °C
Length x Diameter (mm)	106 x 12	140 x 12	140 x 12	114 x 13	140 x 13	140 x 12
Note:	Pin connector, remote fill with side arm, use with indicating half cell	ORP and titrations	Silver and halide titrations	Chlorine titrations with Cl titrimer. Connect to BNC using 13-620-488 (page 35)	KF, dead-stop titrations, sulfur	BNC, most ORP measurements, includes bulb protector

Fisher Scientific accumet Ion-Selective Electrodes (ISE)

Combination And Half Cell ISE's Have BNC For Universal Meter Use

Polymer Membrane And Solid State ISE

Catalog No.	13-620-536	13-620-534	13-620-532	13-620-521	13-620-525
Parameter	Calcium	Nitrate	Potassium	Bromide	Bromide
Type	ISE - polymer membrane	ISE - polymer membrane	ISE - polymer membrane	ISE - solid-state	ISE - solid-state
Combination Or Half Cell	Combination	Combination	Combination	Sensing half cell	Combination
Refillable Or Gel (Sealed)	Refillable	Refillable	Refillable	n/a	Refillable
Refill Solution	13-642-550	ACCU0807-500	Fill solution (Dilute ISA from 5M to 0.1M)	n/a	ACCU0834-500
Ionic Strength Adjuster	13-641-851	13-641-850	13-641-927	ACCU0820-500	13641851
Glass Or Plastic Body	Plastic	Plastic	Plastic	Plastic	Glass
Max Temp	40 °C	40 °C	40 °C	80 °C	80 °C
Length x Diameter (mm)	102 x 13	102 x 13	102 x 13	102 x 13	108 x 13
Connection	BNC	BNC	BNC	BNC	BNC
Interferences	Mg ⁺⁺ , Zn ⁺⁺ , Ba ⁺⁺ , K ⁺ , Na ⁺ , Ni ⁺⁺ , Cu ⁺⁺ , Fe ⁺⁺ , Sr ⁺⁺ , H ⁺⁺ , Hg ⁺⁺ , Pb ^v	Cl ⁻ , NO ₂ ⁻ , Br ⁻ , CN ⁻ , ClO ₃ ⁻ , I ⁻ , ClO ₄ ⁻	Cs ⁺ , NH ₄ ⁺ , Tl ⁺ , H ⁺ , Ag ⁺ , Tris, Li ⁺ , Na ⁺	I ⁻ , CN ⁻ , S ⁻	S ⁻ , I ⁻ , CN ⁻ , High Cl ⁻ levels, NH ₃
Range (ppm)	0.2 to 40,000	0.5 to 14,000	0.04 to 39,000	0.4 to 79,000	0.4 to 79,000
Calibration Solutions	100 ppm 13-641-862	1000 ppm = 13-641-910 100 ppm = 13-641-924	1000 ppm = 650016 100 ppm = 649716	1000 ppm ACCU0822-500	1000 ppm ACCU0822-500
0.1 M	CaCl ₂ (13-641-811)	NaNO ₃ (13-641-888)	KCl (13-641-917)	NaBr (13-641-866)	NaBr (13-641-866)
Typical Applications & Notes:	Foods, beverages, soil, pharmaceuticals, explosives, fertilizers, plants, EDTA titration endpoint	Pollution testing, foods, pharmaceuticals, fertilizers, plants, meats, pickling baths	Body fluids, soils, sewage, fertilizers, foods, beverages	Biological fluids, soil, plants, foods, effluents, Method ASTM approved. Requires reference half cell	Biological fluids, soil, plants, foods, effluents, Method ASTM approved

Solid State ISE

Catalog No.	13-620-627	13-620-519	13-620-547	13-620-629	13-620-523
Parameter	Chloride	Chloride	Cupric	Fluoride	Fluoride
Type	ISE - solid-state	ISE - solid-state	ISE - solid-state	ISE - solid-state	ISE - solid-state
Combination Or Half Cell	Combination	Sensing half cell	Combination	Combination	Sensing half cell
Refillable Or Gel (Sealed)	Refillable	Sealed	Refillable	Refillable	Sealed
Refill Solution	13-620-432	n/a	ACCU0834-500	13-620-431	n/a
Ionic Strength Adjuster	13-641-852	13-641-852	13-641-852	TISAB II 13-642-578 TISAB III 13-641-874	TISAB II 13-642-578 TISAB III 13-641-874
Glass Or Plastic Body	Plastic	Plastic	Glass	Plastic	Plastic
Max Temp	80 °C	80 °C	80 °C	80 °C	80 °C
Length x Diameter (mm)	110 x 13	102 x 13	108 x 13	110 x 13	102 x 13
Connection	BNC	BNC	BNC	BNC	BNC
Interferences	S ⁻ , I ⁻ , CN ⁻ , OH ⁻ , Br ⁻	Br ⁻ , I ⁻ , CN ⁻ , S ⁻ , OH ⁻	Ag ⁺ , Hg ⁺⁺ , Cl ⁻ , Br ⁻ , Fe ⁺ , Cd ⁺⁺	OH ⁻	OH ⁻
Range (ppm)	1.8 to 35,500	1.8 to 35,500	0.00064 to 6,350	0.02 to saturated	0.02 to saturated
Calibration Solutions	1000 ppm (13-299-87)	1000 ppm (13-299-87)	1000 ppm (23004)	100 ppm (13-641-872) 1 ppm w/TISAB II (15-411) 2 ppm w/TISAB II (15-412) 10 ppm w/TISAB II (13-641-752)	100 ppm (13-641-872) 1 ppm w/TISAB II (15-411) 2 ppm w/TISAB II (15-412) 10 ppm w/TISAB II (13-641-752)
0.1 M	NaCl (13-641-876)	NaCl (13-641-876)	Cu(NO ₃) ₂ (13-641-835)	NaF (13-641-871)	NaF (13-641-871)
Typical Applications & Notes:	Water/wastewater, soil, dairy, meats, tomato/vegetable products, Method ASTM/AOAC approved	Water/wastewater, soil, dairy, meats, tomato/vegetable products, Method ASTM/AOAC approved. Requires reference half cell such as 13-620-216.	Plating baths, water	Drinking water, stack gases, sea water, minerals, soils, foods, biological fluids, toothpaste, ASTM/EPA Method	Drinking water, air and stack gases, sea water, minerals, soils, foods, biological fluids, toothpaste

Solid State ISE

Catalog No.	13-620-549	13-620-543	13-620-551	13-620-545	13-620-538
Parameter	Lead	Lead	Silver/sulfide	Silver/sulfide	Cyanide
Type	ISE - solid-state	ISE - solid-state	ISE - solid-state	ISE - solid-state	ISE - solid-state
Combination Or Half Cell	Combination	Sensing half cell	Combination	Sensing half cell	Combination
Refillable Or Gel (Sealed)	Refillable	n/a	Refillable	n/a	Refillable
Refill Solution	ACCU0834-500	n/a	ACCU0834-500	n/a	ACCU0808-500
Ionic Strength Adjuster	-	-	13-641-852 SAOB 13-641-882	13-641-852 SAOB 13-641-882	13-641-853
Glass Or Plastic Body	Glass	Plastic	Glass	Plastic	Plastic
Max Temp	80 °C	80 °C	80 °C	80 °C	80 °C
Length x Diameter (mm)	108 x 13	102 x 13	108 x 13	108 x 13	102 x 13
Connection	BNC	BNC	BNC	BNC	BNC
Interferences	Ag ⁺ , Hg ²⁺ , Cu ²⁺ , CD ²⁺ , Fe ²⁺	Ag ⁺ , Hg ²⁺ , Cu ²⁺ , CD ²⁺ , Fe ²⁺	Hg ²⁺	Hg ²⁺	Cl ⁻ , Br ⁻ , I ⁻ , S ²⁻ absent
Range (ppm)	0.2 to 20,700	0.2 to 20,700	0.01 to 107,900 (Ag ⁺) 0.003 to 32,100 (S ²⁻)	0.01 to 107,900 (Ag ⁺) 0.003 to 32,100 (S ²⁻)	0.1 to 260
Calibration Solutions	100 ppm (13-641-774) Pb(ClO ₄) ₂ (13-641-773)	- Pb(ClO ₄) ₂ (13-641-773)	-	-	-
0.1 M	-	-	-	-	-
Typical Applications & Notes:	Organic compounds, water/wastewater, plating baths	Organic compounds, water/wastewater, plating baths. Requires reference half cell	Sewage effluent, soils, sediments, plating baths, pulping liquors, photographic fixing solution	Sewage effluent, soils, sediments, plating baths, pulping liquors, photographic fixing solution. Requires reference half cell	Silicon petrochemical, plating water, wastes

Gas-Sensing And Glass Membrane ISE

Catalog No.	13-620-509	13-620-511	13-620-503A	13-620-501
Parameter	Ammonia	Carbon Dioxide	Sodium	Sodium
Type	ISE - gas sensing	ISE - gas sensing	ISE - glass membrane	ISE - glass membrane
Combination Or Half Cell	Combination	Combination	Combination	Sensing half cell
Refillable Or Gel (Sealed)	Refillable	Refillable	Refillable	Refillable
Refill Solution	ACCU0800-500	ACCU0806-500	SP138-500	SP138-500
Ionic Strength Adjuster	13-641-883	13-641-770	13-641-748	13-641-748
Glass Or Plastic Body	Plastic	Plastic	Plastic	Glass
Max Temp	50 °C	50 °C	60 °C	80 °C
Length x Diameter (mm)	108 x 12	108 x 12	106 x 12	108 x 13
Connection	BNC	BNC	BNC	BNC
Interferences	Volatile amines, metal cations that complex ammonia	Volatile organic acids	Ag ⁺ , Li ⁺ , K ⁺ , NH ₄ ⁺	Ag ⁺ , Li ⁺ , K ⁺ , NH ₄ ⁺
Range (ppm)	0.009 to 1,700	0.44 to 1,320	0.023 to 23,000	0.023 to 23,000
Calibration Solutions	100 ppm (13-299-98) 1000 ppm (13-641-924C)	1000 ppm (13-641-769)	10 ppm (13-641-948) 100 ppm (13-641-909) 1000 ppm (13-641-747) 1000 ppm w/ISA (13-641-787)	10 ppm (13-641-948) 100 ppm (13-641-909) 1000 ppm (13-641-747) 1000 ppm w/ISA (13-641-787)
0.1 M	NH ₄ Cl (13-641-923)	NaHCO ₃ (13-641-768)	10 % NaCl (13-641-876)	10 % NaCl (13-641-876)
Typical Applications & Notes:	Sewage effluent, boiler water, fish tanks, stack gases, sea water, biological samples, fertilizers. Uses 13-620-512 replacement membranes	Measures carbon dioxide, carbonate, bicarbonate in soft drinks, wines, ground/sea water, fermentation. Uses 13-620-513 replacement membranes	Steam condensate, meats, fish, dairy, fruit juices, fermentation, ground/sea water, soils, body fluids	Steam condensate, meats, fish, dairy, fruit juices, fermentation, ground/sea water, soils, body fluids. Requires reference half cell



Fisher Scientific accumet Conductivity, Dissolved Oxygen And Temperature Probes

Keep Your Instruments Running Smoothly

Conductivity Cells

Conductivity/Temperature Electrodes For Fisher Scientific accumet™ AB200, XL200, XL500, And XL600 Benchtop Meters

Sensor Type	Body Type	Nominal Cell Constant	Ideal Range	Catalog No.
2-Cell	Epoxy	0.1	0.5 to 2,000 μ S	13-620-101
2-Cell	Epoxy	1	20 μ S to 20 mS	13-620-100
2-Cell	Epoxy	10	1 to 200 mS	13-620-102
4-Cell	Glass	1	10 μ S to 100 mS	13-620-163
4-Cell	Glass	10	1 to 200 mS	13-620-164
4-Cell	Epoxy	1	10 μ S to 100 mS	13-620-165
4-Cell	Epoxy	10	1 to 200 mS	13-620-166

Cell For Fisher Scientific accumet™ AP75 And AP85 Portable Meters

2-Cell	Ultem	1	20 μ S to 20 mS	13-620-AP54
--------	-------	---	---------------------	--------------------



Dissolved Oxygen Probes

Dissolved Oxygen Probes And Accessories For Fisher Scientific accumet™ AR60, AR40, AB40, XL40, XL60, And XL600 Benchtop Meters

Description	Catalog No.
Self-Stirring DO/BOD/Temp Probe	13-620-SSP
Adapter To Connect YSI™ Self Stirring DO Probes To AR, AB, XL Series Meters	13-637-DOADPT
Membrane Kit For 13-620-SSP. Includes (6) Membrane Caps, Polishing Disk, And Electrolyte Filling Solution	13-637-DOM

Dissolved Oxygen Probes And Accessories For Fisher Scientific accumet™ AP74 And AP84 Portable Meters

Description	Catalog No.
DO/Temperature Probe With 10-Ft Submersible Cable	15-500-034
DO/Temperature Probe With 25-Ft Submersible Cable	15-500-035
Electrolyte Solution For Dissolved Oxygen Probes, 500 mL	15-500-036
DO Replacement Membrane Assembly, Preassembled Membrane, Membrane Lock, O-ring, Cap	15-500-037
Replacement DO Probe Membranes, Pack Of 25, Requires Membrane Installation Tool	15-500-038
Membrane Installation Tool, Required To Replace Membranes	15-500-039



ATC/Temperature Probes

All temperature probes are stainless steel and have a 3-ft cable.


Description	Catalog No.
ATC For XL, AR, And AB Series Benchtop Meters	13-620-19
ATC For AP72 And AP71 Portable Meters	13-620-20
ATC For AP100 And AP60 Series Portable Meters	13-620-AP53
ATC For Discontinued Fisher Scientific accumet Models 10, 15, 20, 25, 30, 50, 150, 750, 800MP Series, And 900 Series	13-620-16





BNC Adapter Cables

Description	Catalog No.
U.S. Standard/Pin To BNC Input Jack	13-620-498
Pin Plug To BNC Input Jack. Use With Metallic Electrodes Having Pin Connectors	13-620-490
BNC/Pin Plugs To Single BNC Input Jack	13-620-489
Dual-Pin Plugs To BNC Input Jack	13-620-488

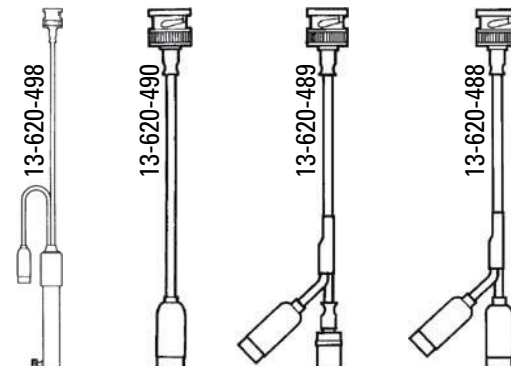

Precision Calibration Resistor Kit

Check Meter Compliance To USP Standards. Includes:

- Six Precision Resistors (Nominal Values 10 ohm, 100 ohm, 1 kilohm, 10 kilohm, 100 kilohm, And 1 megohm)
- Connectors For Fisher Scientific accumetXL, AR, And AB Conductivity Meters
- Foam Padded Case

Description	Catalog No.
Fisher Scientific accumet Conductivity Calibration Kit	13-637-674



Need this NIST certified? Ask for Innocal 17090-30 certificate on [page 47](#).



Other Replacement Parts And Accessories

Description	Catalog No.
Fisher Scientific accumetR Benchtop Stirring Probe	13-620-BSP
Replacement Paddle For 13-620-BSP	13-620-RP
Compact Thermal Printer For ABxxx And XLxxx Series, 100-240 V. Includes (2) Rolls Paper And 4 Plug Types For Global Use	13-637-690
Replacement 5-Yr Thermal Paper For 13-637-690 (Pack Of 2 Rolls)	13-637-691
Electrode Arm And Bracket For All AB And XL Series Benchtop Meters	13-637-671
Fisher Scientific Optical USB Mouse. Compatible With Any XL Series	13-637-692
Replacement Power Supply For AB150, AB200, AB250, XL150, XL200, XL250, XL500, XL600, 100/240 VAC	13-636-104
Replacement RS-232 Output Cable For PC And Printer	13-637-681



Certified Buffer Solutions

Standardized Against NIST Standard Reference Material

Thermo Scientific™ Orion™ Certified Color-Coded pH Buffers



- Tight tolerances:
 - 0.01 for pH 4.0 and 7.0
 - 0.02 for pH 10.0
- Standardized at 25 °C against NIST standard reference material
- Label includes temperature correction chart

pH	Color Coded	Size	Catalog No.
1.68		475 mL	13-643-002
4.01	Red	475 mL	13-301-133
5.00	Orange	475 mL	13-640-8
6.86		475 mL	13-643-003
7.00	Yellow	475 mL	13-641-857
9.18		475 mL	13-643-004
10.01	Blue	475 mL	13-079-319
12.46		475 mL	13-643-001

For best results, always use fresh buffer solutions. Don't use expired buffers!



Oakton™ Conductivity Solutions

- Choose 1 pint (475 mL) plastic bottles or (20) single-use pouches

Value	Quantity	Catalog No.
10 µS	20 pouches	50-632-479
23 µS	475 mL	13-300-114
84 µS	475 mL	13-300-111
447 µS	475 mL	13-300-117
447 µS	20 pouches	13-300-129
1413 µS	475 mL	13-300-112
1413 µS	20 pouches	13-300-130
1500 µS	475 mL	13-300-110
2070 µS	475 mL	13-300-115
2764 µS	475 mL	13-300-113
2764 µS	20 pouches	13-300-131
8974 µS	475 mL	13-300-119
12880 µS	475 mL	13-300-109
15000 µS	475 mL	13-300-118
15000 µS	20 pouches	13-300-132
80000 µS	475 mL	13-300-116



pH Buffers For Low Ionic Strength Samples. Compatible With XL Pure Water Buffer Group



13-641-164

13-641-165

13-641-722

Contact our team of knowledgeable electrochemistry experts for recommendations, troubleshooting, and more!
888-358-4706 or
accumet@fishersci.com



Description	Catalog No.
Pure Water Low Ionic Strength pH Test Kit – 4 x 475 mL of pH 6.97 Buffer, 2 x 475 mL Of pH 4.10 Buffer, 2 x 60 mL Of pH-ISA Adjuster	13-641-735
Pure Water pH-ISA Low Ionic Strength Adjuster, 5 x 60 mL	13-641-722
Pure Water pH 4.10 Buffer For Low Ionic Strength pH Samples, 4 x 475 mL	13-641-165
Pure Water pH 6.97 Buffer For Low Ionic Strength pH Samples, 4 x 475 mL	13-641-164
Pure Water pH 9.15 Buffer For Low Ionic Strength pH Samples, 4 x 475 mL	13-641-166

Other Solutions and Accessories

Description	Catalog No.
Saturated KCl electrode filling solution. Use With Double-Junction And Calomel pH Electrodes (Purple Ring), Or Fluoride (13-620-529), Sodium, And Calcium Combination Ion Selective Electrodes, 500 mL	SP138-500
4 M KCl Saturated with AgCl electrode filling solution. Use with Single-Junction pH Electrodes (Blue Ring), 500 mL	SP135-500
Skylite Electrolyte with Syringe for accuCap 13-620-130 pH Electrode, 60 mL	13-636-430
pH Electrode Storage Solution, 500 mL	13-300-178
Electrode Storage Bottle	13-620-499
Deionized Water Rinse Solution, Box of (20) single-use pouches	13-300-180
ORP Standard, +220 mV vs. Ag/AgCl electrode using 13-641-686 fill solution, 475 mL	13-641-210
Electrode storage sleeve and base	13-642-145

13-620-499



13-641-210



13-636-430





About pH, ORP, And ISE Measurement

A Practical Approach To Achieve Best Results


Next to temperature, pH is quite possibly the most common laboratory measurement today. It crosses over many disciplines from water/wastewater, R&D, environmental, chemical & life sciences, and an endless number of industrial applications.

The pH or **p**ower of **H**ydrogen, is the degree of acidity or alkalinity of a solution based on the hydrogen ion activity, represented by the equation:

$$\text{pH} = -\log [\text{H}^+]$$

Stated mathematically, pH is the negative logarithmic value of the Hydrogen ion. As it is based on a log scale, each pH unit represents a factor of 10, so a solution with a pH of 5 is 100 times more acidic than pH of 7.

Here are examples of pH in a few common industrial and household products.

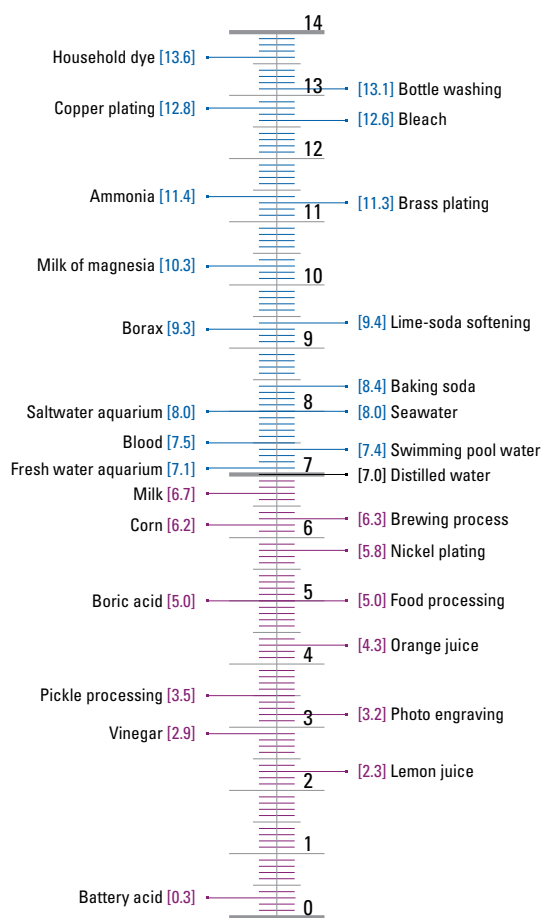

Litmus paper can be used for rough pH measurements, but for highest accuracy a potentiometric system using a pH meter and pH electrode is used. The electrode is sensitive to H⁺ ions, which create a small voltage potential, which is converted to a pH value by the pH meter. Electrode behavior is described by the Nernst equation:

$$E = E_o + (2.3 \text{ RT/nF}) \log a_{\text{H}^+}$$

E is the measured potential from the sensing electrode, E_o is related to the potential of the reference electrode, (2.3 RT/nF) is the Nernst factor and log a_{H⁺} is the pH. The Nernst factor, 2.3 RT/nF, includes the Gas Law constant (R), Faraday's constant (F), the temperature in degrees Kelvin (T) and the charge

of the ion (n). For pH, where n = 1, the Nernst factor is 2.3 RT/F. Since R and F are constants, the factor and therefore electrode behavior is dependent on temperature.

The pH electrode system consists of two half cells: a pH indicating electrode, which develops a potential dependent on the pH of a solution; and a reference electrode, which provides a constant potential and completes the electrical circuit. Using separate pH indicating and reference half cells allows you to select each cell without compromise, tailor the system precisely to the sample's nature, and achieve good accuracy. It can mean lower replacement costs too, since usually only one of a pair is broken. For ATC, a third electrode is required. Nevertheless, the combination electrode – an indicating half cell and a reference half cell joined co-axially – is by far the most common choice for the convenience and compactness it offers. Some combined electrodes also offer built-in ATC for added convenience.

» Temperature Compensation

In a perfect pH electrode – one that measures zero mV at exactly pH 7 – there is no temperature effect on the electrode sensitivity at pH 7 regardless of temperature change. While pH electrodes are not perfect, errors related to temperature are negligible near pH 7, and can be disregarded. As a rule however, the further from pH 7 the solution is and the greater the temperature changes, the greater the expected measurement error due to changes in the electrode's sensitivity. For most electrodes, the error is approximately 0.003 pH/°C/pH away from pH 7.

Consider a pH meter that was calibrated at room temperature (25 °C) and measures a sample at pH 4 at 5 °C,

Temperature difference : 25 °C - 5 °C = 20 °C
 Number of pH units away from neutral : 7 pH - 4 pH = 3 pH units
 Total error : 0.003 x 20 x 3 = 0.18 pH

pH meters that incorporate Automatic Temperature Compensation (ATC) are able to overcome this error.

Q: So why does my pH 10.00 buffer read as 10.06?

A: ATC or not, temperature influences the pH of all solutions. While pH 10.00 buffer is certified to 10.00 at 25 °C / 77 °F, at 20 °C / 68 °F which is often room temperature, it is actually 10.06.

To understand the pH/temperature relationship, we need to look at two main processes; Calibration & Measurement.

Calibration: Meters have a buffer/temperature look-up table stored into memory for every pH buffer that the meter can calibrate to. For example, when calibrating with a pH 10 buffer that is 20 °C, the meter knows to calibrate to a value of 10.06 when calibration is performed. Had the temperature of the buffer been 25 °C, the meter would calibrate instead to 10.00. Using a temperature probe during calibration allows best accuracy as the meter is able to calibrate to the most appropriate pH value.

Measurement: As temperature changes, the pH electrode responds slightly differently. Since this change is predictable, the meter is able to make the appropriate adjustment. The meter senses the temperature and applies the appropriate electrode slope (mV per pH unit) for that temperature. For example, if the temp is 25 °C, the instrument uses 59.16 mV for each pH unit. If the temperature is 20 °C, the meter uses 58.16 mV for each pH unit. Using a temperature probe during measurement allows best accuracy as the meter is able to correct for changes in electrode response due to temperature. Note: Temperature compensation only corrects for the change in the output of the electrode, not for the change in the actual solution pH. As such, the AB15 will always display the actual pH of the solution at the current temperature.

» pH Standardization Buffers

Buffers – solutions of known pH value – allow adjustments to the meter/electrode system to reflect accurate measurements. Certified accurate buffers are available as ready-to-use color coded solutions, concentrated solutions, capsules and prepackaged salts. All have the special characteristic of resisting pH change upon dilution or acid/base contamination. For best accuracy, use a minimum of two-point standardization; first with a buffer value close to the electrode system's zero potential (typically pH 7); and next with an acid or base buffer whose value brackets the expected pH value of the sample. Microprocessor-based meters may permit additional calibrations – up to five points in some models. For best accuracy, perform calibration with ATC at the same temperature as the expected samples.

» ORP Measurement

Oxidation-Reduction Potential (ORP) or “Redox” measurements are used to monitor chemical reactions, to quantify ion activity, or to determine the oxidizing or reducing properties of a solution. ORP is a measurement of the total electrical potential of a solution – a sum of all oxidation and reduction reactions.


ORP electrodes measure the voltage across a circuit formed by the measuring metal half cell and the reference half cell. When the ORP electrode is placed in the presence of oxidizing or reducing agents, electrons are constantly transferred back and forth on its measuring surface, generating a tiny voltage. ORP is expressed as +/- mV, and can be measured by using the millivolt mode of any pH meter with an ORP electrode.

Like pH, an ORP electrode system consists of a sensing and reference electrodes. For ORP, the indicating electrode is metallic – usually platinum, gold, or silver.

ORP varies with temperature, but unlike pH, it is not completely predictable and therefore, ATC is not used. Constant temperature is best for monitoring ORP.

In principle, ORP measurements should not require standardization; in practice however, it may be necessary to check the system against standards of known potential, as described in ASTM Method D 1498.

ORP is also useful in pool water treatment as an indication of sanitation in relation to free chlorine. ORP technology has gained recognition worldwide and is found to be a reliable indicator of bacteriological water quality. The table below illustrates the Kill Time of E.Coli bacteria as a function of ORP value. With a value of 600 mV, the life of the bacteria is almost 2 minutes; at 650 mV it reduces to 30 seconds. Above 700 mV the bacteria is killed within a few seconds. It is therefore necessary for the water to have an ORP value of at least 700 mV to ensure good water quality.

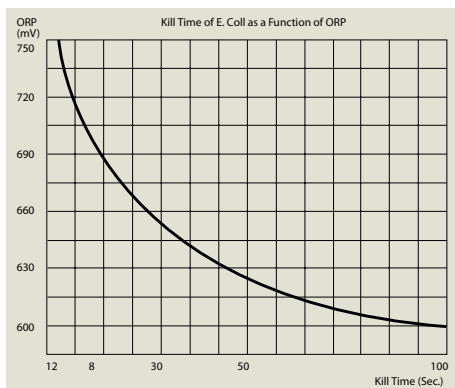

ORP value also depends on the pH of pool water. The pH of the pool water has to be maintained at the optimum level between 7.2 and 7.6 pH by dosing appropriate chemicals. If the pH of the swimming pool water is acceptable and the ORP value is below 700 mV, hypochlorite or other oxidizing chemicals should be added.

Common ORP applications include the treatment of industrial wastes, study of biological systems, oxidation of cyanide, bleaching of pulp, manufacture of bleach and reduction of chromate wastes.

» ISE Measurement

Ion-Selective Electrodes (ISE) respond to a particular ion in solution because of the characteristics of the electrode's sensing membrane. Ideally, the ISE develops an electrical potential which is proportional to the concentration of the ion for which the membrane is selective. The most widely-used ISE is the glass-membrane pH electrode (selective for hydrogen ion).

When an ISE – the indicator electrode – and a reference electrode are placed in a solution and connected to a pH/ion meter, they form a potentiometric cell. At equilibrium, the meter measures the potential difference between the ISE and the reference electrode. This potential is proportional to the activity of the ion of interest.

The table below gives theoretical slope values at 25 °C:

Species	Slope (mV/Decade)
Monovalent Cation	+59.16
Monovalent Anion	-59.16
Divalent Cation	+29.58
Divalent Anion	-29.58

Although ISE's are most often measured by direct analysis, some sample preparation is required. Normally, an ionic strength adjuster must be added equally to samples and standards. Then, standards must be prepared which are used to calibrate the meter, or to construct a calibration curve (by plotting the electrode's output in mV vs. the log of concentration). Sample concentration can then be read directly from the meter or calibration curve.

Incremental methods can reduce errors caused by temperature variations, complex matrices, and complexation. They're also useful for applications where only occasional samples are analyzed. Incremental methods include:

- 1) Known Addition;
- 2) Known Subtraction;
- 3) Analate Addition; and
- 4) Analate Subtraction.

Ion-selective electrodes can also be used to detect the endpoint of a titration. The ISE can be selected to monitor either the addition of titrant or the depletion of analate. The electrode potential is plotted vs. the volume of titrant added. The volume corresponding to the equivalence point is determined from the graph, and used to calculate sample concentration.

A number of metallic electrodes are also used in titrations. Dual platinum wire and plate electrodes, are used with pH meters and titration instruments in dead-stop and amperometric titrations; and silver billet electrodes are the choice for silver and halide titrations.

The ion meter/electrode system must be standardized by immersing the electrodes into solutions having a known concentration of the ion of interest. Stock ISE standards are available in a variety of molar, ppm, and percent concentrations. Calibration is typically done using at least two points that are 10 fold apart, such as 1 and 10 ppm. An ionic strength adjuster is required to eliminate interference from other ions. Since some ISE's have a restricted pH range, a pH adjustment solution may also be necessary. Other special reagents and solutions are available for specific applications.

Testing/Care/Troubleshooting


Prevent Problems From Happening And Recognize When It's Too Late

» pH Electrode Offset

A perfect pH electrode under ideal conditions, when placed in pH 7.00 buffer at 25 °C, will produce 0 mV. The difference between 0 mV and the electrode's actual mV reading is called the offset error, which usually becomes problematic when it is more than ±30 mV or about ½ of a pH unit.

» pH Electrode Slope

A perfect pH electrode under ideal conditions at 25 °C, produces a slope of 59.16 mV per pH unit. For example, an electrode with 0 mV offset should read mV value of 177.48 mV when placed in a pH 4.01 solution. The slope is calculated as $(177.48 \text{ mV} - 0 \text{ mV}) / 3 \text{ pH} = 59.16 \text{ mV/pH}$. The difference between this perfect slope reading and the electrode's actual reading is called the slope error. These theoretical values are not always achieved, even with brand new electrodes. The slope of most new pH electrodes should fall between 92 % and 102 % of 59.16 mV.

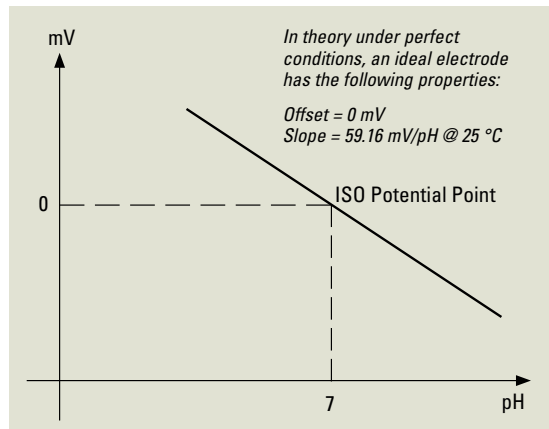

» Normal Aging

As electrodes are used or stored for long periods they can undergo changes in performance. The electrode offset and slope can be monitored to evaluate these changes. Periodic calibration is usually adequate to correct for these changes. If an electrode is able to be calibrated and reads accurately in certified standards, is stable, is responsive and repeatable, it is still considered functional and may be used indefinitely although it no longer meets "new" electrode specifications with regards to offset or slope.

» How to check your pH electrode – including slope and offset:

1. Clear/reset any existing calibration stored in your meter then select the millivolt (mV) mode of your pH meter
2. Using a pH electrode, obtain mV readings of two fresh calibration buffers (i.e. pH 4.0 and pH 7.0 are best)
3. Determine the net mV change
4. Determine the net mV change per pH unit change and compare using the chart below. For example, at 25 °C: pH 4 = 170.5 mV, pH 7 = -3.4 mV: Net change = $170.5 \text{ mV} - (-3.4 \text{ mV}) = 173.9 \text{ mV}$. Since pH 4 and pH 7 are 3 pH units apart, using the chart below, 98 % = Very Good
5. The offset can be determined by observing the mV reading with your pH electrode in pH 7 buffer. The closer to 0 mV the better. A mV more than 30 mV from 0 may be problematic.

6. Also take notice of the electrode response time – the faster, the better. Don't expect an economy gel electrode to respond as fast as a high performance glass refillable electrode, but note that a great slope alone is not meaningful if the electrode takes 2 hours to stabilize.

Slope	At 20 °C		At 25 °C		Status
	1 pH Unit	3 pH Units	1 pH Unit	3 pH Units	
>102	>59.3	>178.0	>60.3	>181.0	Poor
102 %	59.3	178.0	60.3	181.0	Ok
101 %	58.8	176.3	59.8	179.3	Very Good
100 %	58.2	174.5	59.2	177.5	Ideal
99 %	57.6	172.8	58.6	175.7	Very Good
98 %	57.0	171.0	58.0	173.9	Very Good
97 %	56.4	169.3	57.4	172.2	Very Good
96 %	55.8	167.5	56.8	170.4	Very Good
95 %	55.3	165.8	56.2	168.6	Very Good
94 %	57.0	171.0	58.0	173.9	Ok
93 %	54.1	162.3	55.0	165.1	Ok
92 %	53.5	160.5	54.4	163.3	Ok
<92	<53.5	<160.5	<54.4	<163.3	Poor

» Common Symptoms/Cause (Remedy)

• No Response, All Buffers Or Samples Read The Same pH – Usually pH 7.00 Or 0 mV

Broken sensing bulb or wiring problem (replace electrode), probe not connected to input (verify correct channel selection when using multiple-channel meters), probe is not in contact with sample (remove electrode storage bottle or rubber bulb guard), meter automatically has frozen reading (verify that the hold feature or auto read feature is set to off when using meters with this feature, usually by pressing measure or read.)

• Slow Response With Excessive Crystallization Inside Probe Of Refillable Electrodes

Electrolyte flow clogged from supersaturated electrolyte ("flush & fill" by remove the filling solution through the fill hole with a syringe or by shaking it upside down. Repeatedly flush and rinse the reference cavity with clean, 60-80 °C water to dissolve crystals until removed. Replace filling solution and apply gentle pressure to the filling hole. Re-hydrate electrode in storage solution or pH 4 buffer), (ensure fill hole is in open position). To prevent this in the future, ensure that the re-fill hole is closed when electrode is not in use.

• Slow Response Due To Clogged Junction-Usually With Single-Junction Electrodes

Reaction with silver such as silver sulfide formation or protein deposits which causes a dark spot on the ceramic reference junction (For protein layers prepare a 1 % pepsin solution in 0.1 M of HCl and soak the reference junction for one hour in this solution. Rinse the electrode with distilled water. Alternatively, heat a diluted KCl solution to 60 to 80 °C. Place the sensing part of the electrode into the heated solution for about 10 minutes. Allow the electrode to cool in some unheated KCl solution.

• Dried Salt Deposits Present

Electrolyte residue deposited on electrode surface – often with new electrodes or periods of non-use. (simply dissolve the deposits in warm tap water followed by a brief soak in pH 4 buffer).


- **Slow Response, Noisy, Unstable, Or Erratic Readings**

Sensing bulb dry or dirty (clean electrode with mild detergent & warm water and re-hydrate electrode), temperature may be changing rapidly or electrode may be thermally shocked (allow electrode to reach sample temperature), sample may be non-aqueous (take 30 second readings and soak in pH buffer for one minute between measurements).

For additional support contact your Fisher Scientific sales or customer service representative, call a electrochemistry specialist at 1-888-358-4706, or send an email to: accumet@fishersci.com

» General Rules And Tips

- When using refillable electrodes, open the fill hole during calibration and measurement – but remember to close it afterwards when finished!
- The level of electrolyte in the outer cavity of refillable electrodes should be kept above the level of the solution being measured to prevent reverse electrolyte flow.
- The electrode need only be immersed far enough to cover both the glass pH sensing bulb and reference junction to obtain accurate readings.
- Electrodes perform best when they are hydrated. However, if they dry out they can be reconditioned to normal performance again. Soaking in electrode storage solution helps to optimize and re-establish the thin hydration layer on the sensing bulb that is critical to pH measurement.
- Rinsing the electrode with deionized or distilled water between samples is fine, but storage in deionized or distilled can be detrimental as it will rob critical ions from the sensing bulb. Also, avoid wiping or touching the sensing bulb to maintain the hydration layer and producing any electrical charge.
- Moving or touching the electrode cable may result in unstable readings due to the high impedance (resistance) of the pH glass membrane and introduce noise.
- To eliminate temperature errors associated with the electrode, manual or automatic temperature compensation (ATC) should be used for best accuracy. Since temperature changes pH, the sample temperature should always be noted with pH readings. i.e.) Record results as “pH 8.43 @ 23.2 °C”, instead of “pH 8.43”.
- Always use fresh pH buffers for calibration. Excessive air exposure and sunlight alter the buffers value – especially pH 10.00 buffers which is particularly susceptible to drift.

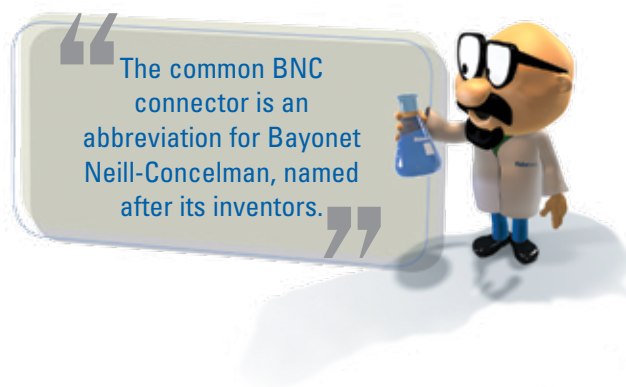

About Conductivity Measurement

A Practical Approach To Achieve Best Results

Electrical Conductivity (EC) meters measure the capacity of ions in an aqueous solution to carry electrical current. As the ranges in aqueous solutions are usually small, the basic units of measurements are milliSiemens/cm (mS) and microSiemens/cm (µS).



Solution	Conductivity
Absolute Pure Water	0.055 µS
Power Plant Boiler Water	1.0 µS
Good City Water	50 µS
Ocean Water	53 mS
Distilled Water	0.5 µS
Deionized Water	0.1 - 10 µS
Drinking Water	0.1 - 1 mS
Wastewater	0.9 - 9 mS
Seawater	53 mS
10 % NaOH	355 mS
10 % H2SO4	432 mS
31 % HNO3	865 mS

» Conductivity Measurement

The principle by which instruments measure conductivity is simple – two plates are placed in the sample, a potential is applied across the plates, and the current is measured. Conductivity (G), the inverse of Resistivity (R) is determined from the voltage and current values according to Ohm's law.


$$G = I/R = I \text{ (amps)} / E \text{ (volts)}$$

Since the charge on ions in solution facilitates the conductance of electrical current, the conductivity of a solution is proportional to its ion concentration. However, conductivity may not correlate directly to concentration. The graphs below illustrate the relationship between conductivity and ion concentration for two common solutions. Notice that the graph is linear for sodium chloride solution, but not for highly concentrated sulfuric acid.

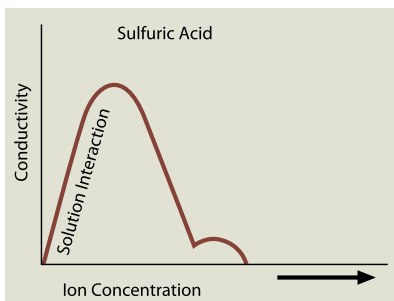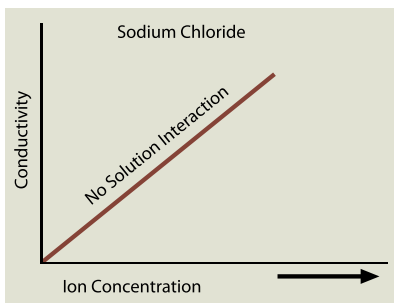

» Resistivity, Conductivity, and Salinity Units

The basic unit of conductivity is the Siemens (S), formerly called the mho. Since cell geometry affects conductivity values, standardized measurements are expressed in specific conductivity units (S/cm) to compensate for variations in electrode dimensions. Specific conductivity (C) is simply the product of measured conductivity (G) and the electrode cell constant (L/A), where L is the length of the column of liquid between the electrode and A is the area of the electrodes.

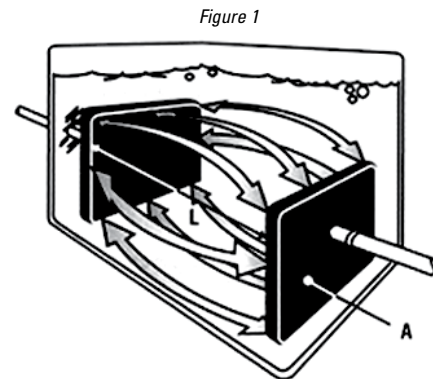

$$C = G \times (L/A)$$

Resistivity is simply the reciprocal of conductivity: conductivity = 1/resistivity. In practice, resistivity units are used when describing ultra pure water such as deionized or reverse-osmosis water. Conductivity units are used for water ranging from drinking water to brackish water. Salinity units are reserved for the highest concentrations expressed as parts per thousand (ppt) or more commonly % concentrations of salt, such as 3.5 % sea water.

Conductivity	Resistivity
0.01 µS	100 MΩ
0.055 µS	18.0 MΩ
0.1 µS	10 MΩ
1 µS	1 MΩ
10 µS	0.1 MΩ
100 µS	0.01 MΩ
1 mS	1 kΩ

» Total Suspended Solids (TSS) and Total Dissolved Solids (TDS)

Suspended Solids measurements are traditionally determined by filtering a measured volume of water – usually 1 liter. Solids captured by a filter are deemed “suspended solids”. The water that passes thru (filtrate) is oven dried and the residue that remains is called “dissolved solids”. Both are usually measured as mg/L or ppm. Unfortunately, drying of the filtrate to complete evaporation is a long process and requires heating. TDS (conductivity multiplied by a TDS factor) can be used as a quick, inexpensive alternative method. TDS as measured with a conductivity meter will correspond to the dissolved ionic content that is often a useful approximation for the total dissolved material.

There are limitations when using TDS. First, a TDS factor used is salt specific so if there are multiple or unknown salts in solution, it's nearly impossible to determine the ideal factor to use. Second, since ionic concentrations are not linear, the TDS factor changes with concentration. Also, TDS is not preferred for low measurements.

» Temperature Compensation and Coefficient

Conductivity is greatly influenced by temperature. Most fluids increase in conductivity as temperature increases. Most ionic solutions will increase about 2 % for each 1 °C increase. Unfortunately, this temperature coefficient (TC) is not always. In the case of high resistance water it can be closer to 5 % or so per °C.

Many instruments adjust the conductivity value based on a TC and display a value that is said to be corrected or normalized to 25 °C. The meter will automatically make corrections to the reading and display a value as if the sample was 25 °C, no matter what the actual temperature is. Some instruments use a fixed TC of 2.0 % per °C. Let us consider a meter that uses 2.0 % TC to measure a 1413 µS standard at 25 °C (77 °F). If the standard is warmed to 30 °C (86 °F), the meter applies a correction of 5 degrees x 0.02 % x 1413 µS = 141.3. Without correction (0.0 % TC) the actual value of a 1413 µS standard of KCl at 30 °C (86 °F) is 1548 µS. As the meter corrects for temperature, it displays a value of 1548 µS minus 141.3 µS = 1407 µS. When the sample cools to 25 °C, it will again read 1413 µS as no correction is applied. Although conductivity cell response is immediate, temperature corrected values will fluctuate as the temperature measurement stabilizes. Advanced meters offer adjustable TC's, usually from 0.0 % to as much as 10 % per °C.


» Calibration and Maintenance

Conductivity systems should be calibrated to certified solution standards before using. When selecting a standard, choose one that has the approximate conductivity of the solution to be measured. Unlike pH, conductivity cells generally change less over time and therefore do not require as frequent calibration.

A polarized or fouled electrode must be cleaned to renew the active surface of the cell. In most situations, hot water with a mild liquid detergent is an effective cleanser. Alcohol easily cleans most organic matter, and chlorine solutions will remove algae, bacteria or molds. To prevent cell damage, abrasives or sharp objects should not be used to clean an electrode. A cotton swab works well for cleaning but care must be taken not to widen the distance of cell.


» 2-Cell Vs. 4-Cell

Most conductivity meters use a 2-cell electrode. The electrode surface is usually platinum, titanium, gold-plated nickel, or graphite. The 4-cell electrode is more expensive and uses a reference voltage to compensate for any polarization or fouling of the electrode plates. The reference voltage ensures that measurements indicate actual conductivity independent of electrode condition, resulting in higher accuracy.

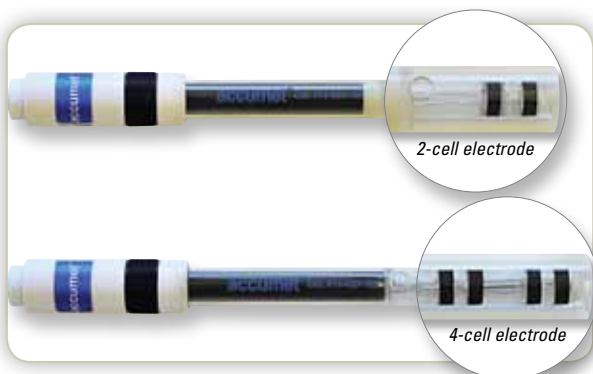

» Advanced Meter Features To Consider

- **Adjustable TDS Conversion Factor**
When a solution does not have a similar ionic content to natural water, the ability to adjust the TDS conversion factor will enable improved accuracy.
- **Adjustable Temperature Coefficients**
Allows for the precise manipulation of the temperature compensation levels for improved results at various temperatures.
- **Multiple Cell Constants**
Allows use of additional electrodes for best results – usually low or high ranges.

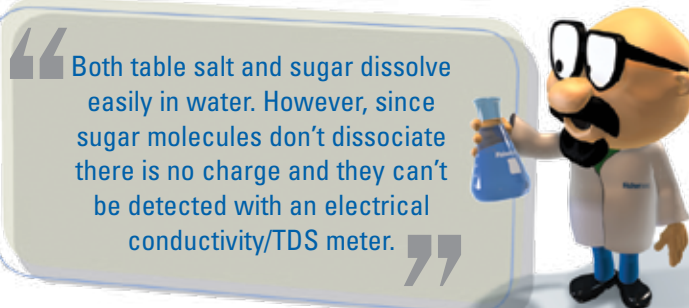

About Dissolved Oxygen Measurement

A Practical Approach To Achieve Best Results

» What is Dissolved Oxygen?

Dissolved Oxygen (DO) is a measure of the amount of dissolved gaseous oxygen in a solution. Some gases, such as ammonia, carbon dioxide and hydrogen chloride, react chemically with water to form new compounds. However, gases such as nitrogen and oxygen merely dissolve in water without chemically reacting with it, and exist as microscopic bubbles between water molecules.

There are two main ways in which dissolved oxygen occurs naturally in water: From the surrounding atmosphere, where oxygen in the surrounding air dissolves readily when mixed into water, up to saturation, during water movements; Via photosynthesis when oxygen is produced by aquatic plants and algae as a by-product of photosynthesis. The amount of oxygen dissolved in water is usually measured in percent saturation, or expressed as a concentration in milligrams per litre water. Accurate measurement of dissolved oxygen is essential in processes where oxygen content affects reaction rates, process efficiency or environmental conditions, such as biological wastewater treatment, wine production, bio-reactions, environmental water testing.

» Basic Principle in DO Measurement

In theory, the amount of DO in a solution is dependent on three factors, namely temperature, salinity and atmospheric pressure.

1. Water Temperature

Solubility of oxygen reduces as temperature increases. Hence, the colder the water, the more dissolved oxygen it contains. Since temperature affects both the solubility and diffusion rate of oxygen, temperature compensation is necessary for any standardized DO measurements.

2. Salinity


The amount of dissolved oxygen increases as salinity level decreases. In other words, freshwater holds more oxygen than saltwater. Since the presence of dissolved salts limits the amount of oxygen that can dissolve in water, the relationship between the partial pressure and concentration of oxygen varies with the salinity of the sample.

3. Atmospheric Pressure

There is a direct proportional relationship between the solubility of dissolved oxygen and the surrounding atmospheric pressure. As pressure decreases with increase in altitude, the amount of dissolved oxygen found in water reduces.

To ensure that your dissolved oxygen is not affected by atmospheric pressure, most instruments include barometric pressure compensation. Advanced meters include a built in barometer that measures and adjusts automatically. However, basic meters will make adjustments only after the user inputs the appropriate pressure manually. Atmospheric pressure correction charts included in with certain instrument manuals are also a helpful reference.

» DO Electrodes

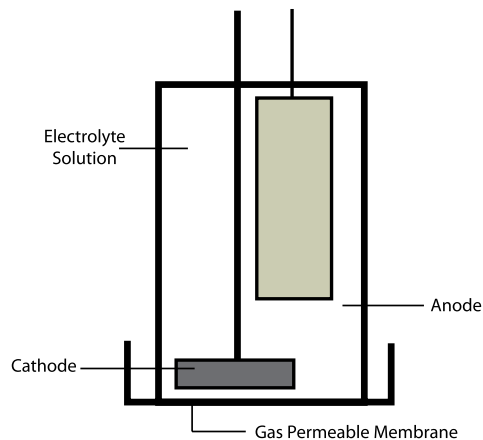

The measurement of DO requires a special DO electrode that is made up of an anode, a cathode, electrolyte solution and a gas permeable membrane. The material of the membrane is specially selected to permit oxygen to pass through. Oxygen is consumed by the cathode which will create a partial pressure across the membrane. Oxygen then diffuses into the electrolyte solution. In short, a DO meter actually measures the pressure caused by movements of oxygen molecules in water or any other medium. The higher the partial pressure of the oxygen in solution, the higher the oxygen concentration. Currently, galvanic and polarographic electrodes are the predominant methods for measuring dissolved oxygen.

The Galvanic Cell consists of two metals, the positive anode and the negative cathode, connected by a salt bridge between the individual half-cells. As the metal electrodes leave electrons behind as they dissolve in the electrolyte. The different properties of the two metals causes them to dissolve at different rates, hence a potential is created when the number of electrons in either side of the cell differs. The potential is translated into an electric current proportion to the oxygen concentration in the electrolyte if an electrical circuit is created between the two electrodes. The galvanic electrode does not need polarising time and is able to assume operation immediately.

During this process, ions of the more active anode are transferred through the electrolyte to the less active cathode, and deposited there as a plating. In this way the anode is corroded. When the anode material eventually corrodes away, the potential drops and the current halts.

The Polarographic Cell consists of two electrodes placed in the electrolyte: One with fixed potential called the reference electrode, and the other with a variable potential called the polarizable electrode. As voltage is applied to the polarizable electrode, a redox reaction occurs, where electrons break away from the electrode to bond with oxygen in the electrolyte. The rate at which the electrons break away from the polarizable electrode is linearly proportionate to the amount of oxygen available in the electrolyte, hence this movement of electrons is representative of the amount of dissolved oxygen left in the electrolyte.

The advantage of a polarographic cell is that the cathode remains intact. The current flow of the polarographic cell is also linearly proportional to the amount of oxygen present in the electrolyte, enabling the cell to provide highly accurate measurements at low oxygen levels.

» BOD & COD

The BOD test measures the molecular oxygen utilized in the biodegradation of organic material and the oxidation of inorganic material. By measuring the amount of oxygen dissolved in samples at the beginning and end of a specified incubation period, the relative oxygen requirements of wastewaters, effluents, and polluted waters can be determined.

$$\text{BOD}_t \text{ (mg/L)} = \frac{D_1 - D_2}{P}$$

BOD_t = Oxygen uptake during incubation period t

D1 = DO of diluted sample immediately after preparation (mg/L)

D2 = DO of diluted sample after incubation period t (mg/L)

P = Decimal volumetric fraction of sample used

BOD is similar to the Chemical Oxygen Demand (COD), which also measures relative oxygen-depletion. However, the possible presence of non-biologically oxidizable may render the COD test to be less accurate.

The COD test is often used to measure the amount of organic compounds in water by measuring the amount of oxygen required to oxidize and break down an organic compound into carbon dioxide, ammonia and water. The basis of the COD test is to determine what can be oxidized into carbon dioxide using a strong oxidizing agent in acidic environments. A blank sample, created by adding all reagents to distilled water is usually used as a control in COD measurements.

Both the BOD and COD tests are means to measure the relative oxygen-depletion effect of a waste contaminant, and are widely used to monitor pollution levels. The BOD test measures the oxygen demand of biodegradable pollutants whereas the COD test measures the oxygen demand of biodegradable pollutants plus the oxygen demand of non-biodegradable oxidizable pollutants.

COD measures everything that can be chemically oxidized and not just the level of biologically active matter that BOD measures. This is especially important to keep in mind when understanding and comparing COD and BOD results that may contain non-biological oxidizable components.






Index

Quick Reference To Every Catalog Number In This Handbook!

Catalog No.	Page	Catalog No.	Page	Catalog No.	Page	Catalog No.	Page	Catalog No.	Page	Catalog No.	Page
13-079-319	36	13-620-164	34	13-620-488	35	13-620-631	15, 26, 28	13-636-AP71A	19	13-641-735	37
13-300-109	36	13-620-165	34	13-620-489	35			13-636-AP72	19	13-641-857	36
13-300-110	36	13-620-166	34	13-620-490	35	13-620-632	25, 26	13-636-AP72A	19	13-642-145	37
13-300-111	36	13-620-181	25	13-620-498	35	13-620-658	30	13-636-AP74	21	13-643-001	36
13-300-112	36	13-620-182A	25	13-620-499	37	13-620-79	30	13-636-AP74A	21	13-643-002	36
13-300-113	36	13-620-183A	15, 25	13-620-501	33	13-620-81	31	13-636-AP75	20	13-643-003	36
13-300-114	36	13-620-185	25	13-620-503A	12, 14, 33	13-620-95	26	13-636-AP75A	20	13-643-004	36
13-300-115	36	13-620-187	25, 29			13-620-96	26	13-636-AP84	21	15-094-26	47
13-300-116	36	13-620-19	15, 34	13-620-509	12, 14, 33	13-620-AP50A	18, 29	13-636-AP84A	21	15-094-27	47
13-300-117	36	13-620-195	24			13-620-AP52	18, 29	13-636-AP85	20	15-094-28	47
13-300-118	36	13-620-196	24	13-620-51	29	13-620-AP53	18, 34	13-636-AP85A	20	15-094-29	47
13-300-119	36	13-620-197	24	13-620-511	33	13-620-AP54	20, 34	13-636-XL150	13	15-094-30	47
13-300-129	36	13-620-20	19, 34	13-620-512	33	13-620-AP55	20, 29	13-636-XL150A	13	15-094-33	47
13-300-130	36	13-620-216	30	13-620-513	33	13-620-AP56	21	13-636-XL200	11	15-094-37	47
13-300-131	36	13-620-221	27	13-620-519	32	13-620-AP61	18, 29	13-636-XL200A	11	15-094-44	47
13-300-132	36	13-620-223A	26	13-620-52	29	13-620-BSP	17, 35	13-636-XL250	12	15-094-50	47
13-300-178	37	13-620-258	30	13-620-521	32	13-620-RP	17, 35	13-636-XL250A	12	15-094-53	47
13-300-180	37	13-620-259	30	13-620-523	32	13-620-SSP	34	13-636-XL250F	12	15-500-034	21, 34
13-301-133	36	13-620-273	30	13-620-525	32	13-621-701	28	13-636-XL250N	12	15-500-035	34
13-620-100	34	13-620-284	29	13-620-53	30	13-621-702	28	13-636-XL500	10	15-500-036	21, 34
13-620-101	34	13-620-285	26	13-620-530A	26	13-621-703	28	13-636-XL500A	10	15-500-037	34
13-620-102	34	13-620-287A	27	13-620-532	32	13-636-100	18	13-636-XL600	9	15-500-038	34
13-620-108A	27	13-620-288	27	13-620-534	12, 14, 32	13-636-104	15, 35	13-636-XL600A	9	15-500-039	34
13-620-111	27, 28	13-620-289	27			13-636-430	37	13-636-XL600D	9	15-700-341	47
13-620-111T	28	13-620-290	27	13-620-536	32	13-636-AB150	15	13-637-671	17, 35	AB150ACERT	47
13-620-112	27	13-620-291	26	13-620-538	33	13-636-AB150A	15	13-637-674	35	AB200ACERT	47
13-620-113	24	13-620-292	26	13-620-543	33	13-636-AB150B	15	13-637-681	35	AB250ACERT	47
13-620-114	24	13-620-293	26	13-620-545	33	13-636-AB200	16	13-637-690	15, 35	SP135-500	37
13-620-115	31	13-620-294	29	13-620-547	32	13-636-AB200A	16	13-637-691	35	SP138-500	37
13-620-122	31	13-620-295	29	13-620-549	33	13-636-AB250	14	13-637-692	35	XL150ACERT	47
13-620-123	31	13-620-296	24	13-620-551	33	13-636-AB250A	14	13-637-DOADPT	34	XL200ACERT	47
13-620-130	25	13-620-297	24	13-620-57	30	13-636-AP110	18	13-637-DOM	34	XL250ACERT	47
13-620-131	25	13-620-298	24	13-620-61	30	13-636-AP110A	18	13-640-8	36	XL500ACERT	47
13-620-132	25	13-620-299A	27	13-620-62	30	13-636-AP115	18	13-641-164	37	XL600ACERT	47
13-620-133	25	13-620-300	27	13-620-627	12, 14, 32	13-636-AP115A	18	13-641-165	37		
13-620-149	31	13-620-31C	28			13-636-AP125	18	13-641-166	37		
13-620-16	34	13-620-45	31	13-620-629	12, 14, 32	13-636-AP125A	18	13-641-210	37		
13-620-163	34	13-620-46	30			13-636-AP71	18	13-641-722	37		

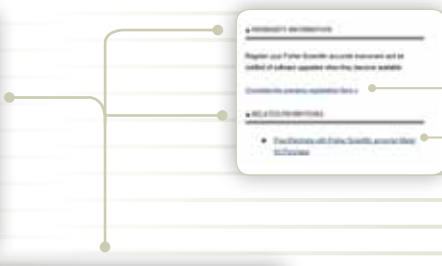

Be sure to visit
www.fishersci.com/accumet
 for product information &
 much more 24 hours a day /
 7 days a week!





INNOCAL[®]
INNOVATIVE CALIBRATION SOLUTIONS

Don't Leave Accuracy To Chance! Order NIST-Traceable Calibration Versions Of Your Favorite Fisher Scientific accumet Meters

Instrument calibration to NIST-traceable standards helps you meet ISO, FDA, USDA, and EPA quality assurance guidelines. Have your Fisher Scientific accumet instrumentation pretested and calibrated before using it in your critical processes.

pH calibration service includes measurement of meter accuracy at five test points across pH range and five test points across corresponding mV range. Testing is performed by our accredited laboratory and an NIST-traceable certificate is provided to document instrument accuracy.



Description	Parameter	Catalog No.	Meter Kit	BioBasic Meter Kit
Fisher Scientific accumet AB Meters				
AB150 with NIST-traceable pH certificate	pH, mV, temp	AB150ACERT	AB150CERT	AB150BCERT
AB250 with NIST-traceable pH certificate	pH, mV, temp, ISE	AB250ACERT	AB250CERT	—
AB200 with NIST-traceable pH certificate	pH, mV, temp, conductivity	AB200ACERT	AB200CERT	—
Fisher Scientific accumet XL Meters				
XL150 with NIST-traceable pH certificate	pH, mV, temp	XL150ACERT	XL150CERT	—
XL200 with NIST-traceable pH certificate	pH, mV, temp, conductivity	XL200ACERT	XL200CERT	—
XL250 with NIST-traceable pH certificate	Dual pH, mV, temp, ISE	XL250ACERT	XL250CERT	—
XL500 with NIST-traceable pH certificate	Dual pH, mV, temp, ISE, conductivity	XL500ACERT	XL500CERT	—
XL600 with NIST-traceable pH certificate	Dual pH, mV, temp, ISE, conductivity, DO	XL600ACERT	XL600CERT	XL600DCERT

NIST-Traceable Certificate For Specifications	Innocal Cert. No.
pH Meter (Benchtop Or Portable) Five test points across pH range and five test points across mV range	17106-20
Conductivity/Resistivity Meter (Benchtop Or Portable) One test point in each measuring range	17090-30
Dissolved oxygen Meter (Benchtop or Portable) Three test points across range	17106-04
Glass Thermometer Two test points across temperature range	17006-03
Humidity Three humidity test points and one temperature test point	17030-20
Infrared Thermometer Four test points across temperature range	17101-62
Pipette/Dispenser Three test points: 25 %, 50 %, and 100 % of range	17104-00
Temperature Datalogger Three test points across temperature range	17002-20
Temperature (Meter Only) Four test points across temperature range	17101-79
Temperature System (Meter + Probe) Four test points across temperature range	17001-10
Timer/Stopwatch Ten test points over a 24-hour period	17060-00



InnoCal is a full-service Metrology Laboratory committed to meeting your calibration needs. Additional instrument testing capabilities are listed below. Contact your Fisher Scientific Customer Service Representative for details on additional calibration capabilities not listed here or call 866-466-6225 to speak with an InnoCal specialist. For calibration service of an instrument you already have, a return authorization number will be issued.

Benefits:

- » **Highest Quality**
Testing performed by an ISO/IEC 17025 accredited laboratory.
- » **Fast Service**
Instruments serviced in 5 to 7 business days.
- » **Excellent Value**
Accuracy verified at multiple test points with "as found" and "as left" test data.
- » **Reliable Documentation**
Certificates are packed with valuable information.

